

## Superior performance in a variety of sterile laboratory filtration applications

- Choose from a variety of membrane types, pore sizes, and effective filtration areas to meet your specific application requirements. Filter sizes accommodate sample volumes from < 2 to 200 mL.
- Improve throughput of viscous or particulate laden solutions with built-in prefilter.

### Product Selection

#### Consider Filtration Volume

The appropriate filter size is determined primarily by the volume being filtered. However, total throughput (capacity) of a filter depends on the characteristics of the fluid being filtered, such as the extent of microbial contamination, viscosity, and particulate load. For small-volume or expensive samples, hold-up volume (sample loss) can be a concern and a smaller filter should be used. The volume recommendations below are intended as a general guide in selecting the most appropriate filter.

Sample Volume (mL)	Filter Size (mm)
up to 10	13
up to 100	25
up to 150	32
up to 200	37

#### Consider Fluid Characteristics

The characteristics of the fluid being filtered are extremely important when evaluating the membrane to be used for filtration. Pall Life Sciences offers sterile Acrodisc syringe filters in the following membrane types:

- Supor® (PES) Membrane**  
For high throughputs, fast filtration, and low protein binding. Not recommended for use with some ketones.
- Versapor® Membrane**  
Economical choice for clarification.
- Fluorodyne® II Membrane**  
For high flow rates and low protein binding. Not recommended for use with some ethers.



- Nylon Membrane**

Used in our DMSO-Safe Acrodisc syringe filter.

- Posidyne® Membrane**

Enhances bioburden and pyrogen removal from aqueous solutions.

- Ultipor® Membrane**

Provides broad solvent and chemical compatibility, and low extractables. Industry standard for pharmaceutical filtration.

#### Consider Scalability

Filters that contain consistent materials of construction provide greatly simplified filter qualification and easier transitions from R&D to pilot to full-scale production. The filters in Pall Corporation's ScaleUp Program incorporate scaleable membranes and materials of construction to assure consistent results, whether you are using syringe filters (less than 100 mL) or larger-capacity capsules and cartridges (up to 5000 L). The ScaleUp Program offers worldwide product availability and validation, troubleshooting, and technical support.

## Sterile Acrodisc Syringe Filters with Supor (PES) Membrane



Universal life science membrane with low protein binding.

- Outstanding flow rate and throughput performance.
- The range of available filter sizes allows you to optimize filtration efficiency without compromising sterilization performance.

### Applications


Use these high-performance filters for prefiltration/clarification of viscous solutions, and sterilization of buffers and culture media. Choose:

- 0.1 and 0.2  $\mu\text{m}$  pore sizes provide sterilization of small volumes of buffers, culture media, and additives.
- Acrodisc PF (0.8 / 0.2  $\mu\text{m}$ ) and Serum Acrodisc syringe filters with unique built-in prefilter for increased throughput of viscous or particulate-laden solutions.
- 0.1  $\mu\text{m}$  pore size filters for reduction of mycoplasma.
- 0.45, 0.8, 1.2 and 5  $\mu\text{m}$  for prefiltration and/or clarification.
- 37 mm Serum Acrodisc GF (Glass Fiber)/0.2  $\mu\text{m}$  for serum and very hard to filter solutions

### Performance


#### Acrodisc and Acrodisc PF Syringe Filters with Supor Membrane

Built-in prefilter enhances throughput of viscous, particulate-laden, or proteinaceous solutions.

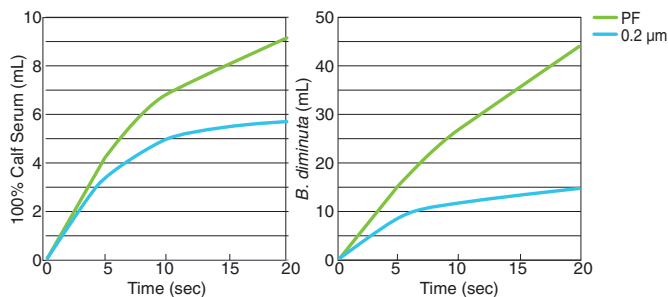


Acrodisc and Acrodisc PF syringe filters with 0.2  $\mu\text{m}$  Supor membrane were challenged with calf serum and a bacterial culture ( $10^7$  cfu/mL) at a constant pressure of 1.4 bar (140 kPa, 20 psi).

## Membrane Filtration Throughput 0.2 $\mu\text{m}$ Rated Membranes

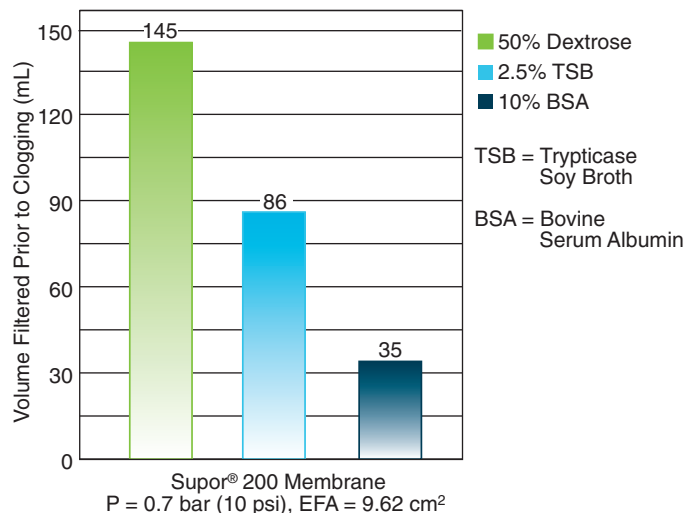

## <sup>125</sup>I BSA Protein Transmission Through 25 mm Acrodisc Syringe Filters

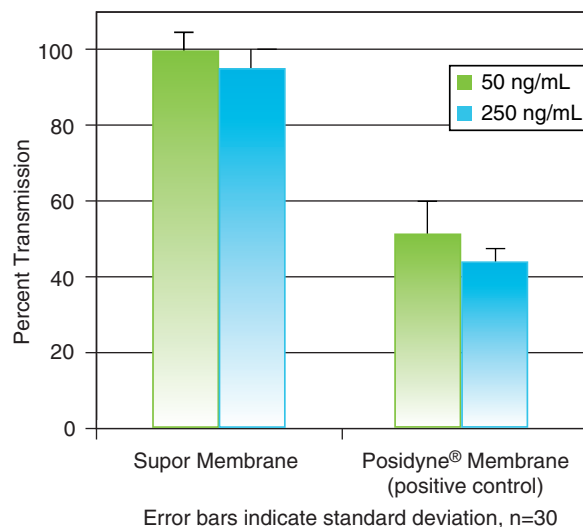

### Specifications

#### Materials of Construction

Filter Media: Supor membrane (hydrophilic polyethersulfone)  
 Serum Acrodisc Prefilter Media: Binder-free borosilicate glass  
 Acrodisc Housing:  
 13 mm: Polypropylene  
 25 and 32 mm: Modified acrylic  
 37 mm: Serum Acrodisc ABS

#### Effective Filtration Area

13 mm: 1.0 cm<sup>2</sup>  
 25 mm: 2.8 cm<sup>2</sup>  
 32 mm: 5.8 cm<sup>2</sup>  
 37 mm: 7.5 cm<sup>2</sup>

#### Inlet/Outlet Connections

Female luer lock inlet, male slip luer outlet

### Typical Hold-up Volume

(with air purge)

13 mm: < 60 µL

25 mm: < 300 µL

32mm: < 500 µL

37mm: < 650 µL

### Maximum Operating Temperature

55 °C (131 °F)

### Maximum Operating Pressure

5.2 bar (520 kPa, 75 psi)

### Typical Water Flow Rate

mL/min at 3.1 bar (310 kPa, 45 psi)

0.1 µm, 25 mm: 35

0.1 µm, 32 mm: > 75

0.2 µm, 13 mm: 22

0.2 µm, 25 mm: 175

0.2 µm, 32 mm: > 330

0.45 µm, 13 mm: 35

0.45 µm, 25 mm: 300

0.45 µm, 32 mm: > 560

0.8 µm, 13 mm: 150

0.8 µm, 25 mm: 700

0.8/0.2 µm, 25 mm: 145

0.8/0.2 µm, 32 mm: > 440

1.2 µm, 32 mm: > 1500

5 µm, 32 mm: > 1725

GF/0.2 µm, 37 mm (Serum Acrodisc Filter): > 425

### Endotoxin

13 & 25 mm : < 0.25 EU/mL using Limulus Amoebocyte

Lysate (LAL) test

32 mm : < 0.4 EU/mL using Limulus Amoebocyte Lysate

(LAL) test

### Sterilization

Sterilized by gamma irradiation and individually packaged

## Ordering Information

### Acrodisc Syringe Filters with Supor Membrane (Sterile)

Part Number	Description	Pkg
4602	0.2 µm, 13 mm	75/pkg
4604	0.45 µm, 13 mm	75/pkg
4608	0.8 µm, 13 mm	75/pkg
4611	0.1 µm, 25 mm	50/pkg
4612	0.2 µm, 25 mm	50/pkg
4614	0.45 µm, 25 mm	50/pkg
4618	0.8 µm, 25 mm	50/pkg
4651	0.1 µm, 32 mm	50/pkg
4652	0.2 µm, 32 mm	50/pkg
4654	0.45 µm, 32 mm	50/pkg
4656	1.2 µm, 32 mm	50/pkg
4650	5 µm, 32 mm	50/pkg

### Serum Acrodisc Syringe Filter (Sterile)

Part Number	Description	Pkg
4525	GF/0.2 µm, 37 mm	20/pkg

### Acrodisc PF Syringe Filters (Sterile)

Part Number	Description	Pkg
4187	0.8/0.2 µm, 25 mm	50/pkg
4658	0.8/0.2 µm, 32 mm	50/pkg

## Acrodisc Syringe Filters Optimised for ScaleUp - Ideal for proof of concept



Four membrane types, Fluorodyne® II, Supor, Posidyne, and Ultipor membranes, assure compatibility with a wide range of fluids.

- Integrity testable (water bubble point).
- Simplify scale up and reduce validation.
- Bacterial retention tested to assure sterile filtrate. Laboratory filters containing the same materials of construction as larger-capacity capsules and cartridges.
- No need to change membrane materials during transition to pilot or full-scale production.

### Applications

- Drug development studies
- Determination of product compatibility and recovery
- Preliminary filterability testing
- Small volume liquid sterilization

### Specifications

#### Materials of Construction

Filter Media: Supor membrane (hydrophilic polyethersulfone), Ultipor membrane (amphoteric Nylon 6,6), Fluorodyne II membrane (hydrophilic polyvinylidene fluoride), Posidyne membrane (positively-charged Nylon 6,6)

Housing: Polypropylene

#### Effective Filtration Area

2.8 cm<sup>2</sup>

#### Inlet/Outlet Connections

Female luer lock inlet, male slip luer outlet

#### Typical Hold-up Volume

(with air purge)

≤ 100 µL

#### Maximum Operating Temperature

60 °C (140 °F) at 2.1 bar (210 kPa, 30 psi)

#### Maximum Operating Pressure

5.4 bar (540 kPa, 80 psi) at ambient temperature

#### Minimum Water Flow Rate

mL/min at 2.1 bar (210 kPa, 30 psi)

PN 4905: 130

PN 4906: 78

PN 4907: 130

PN 4908: 77

#### Recommended Integrity Test

Minimum Bubble Point – Water

PN 4902: 3.32 bar (332 kPa, 48 psi)

PN 4905: 3.5 bar (350 kPa, > 51 psi)

PN 4906: 3.2 bar (320 kPa, > 46 psi)

PN 4907: 3.2 bar (320 kPa, > 46 psi)

PN 4908: 3.2 bar (320 kPa, > 46 psi)

#### Bacterial Retention

Lot samples retain 10<sup>7</sup> cfu/cm<sup>2</sup> of *B. diminuta* per modified

ASTM F838-05

#### Endotoxin

< 0.25 EU/mL using Limulus Amoebocyte Lysate (LAL) test

#### Biological Safety

Passes United States Pharmacopeia (USP) Biological Reactivity Test, *In Vivo* <88>

#### Sterilization

Sterilized by gamma irradiation and individually blister packed

### Ordering Information

#### Acrodisc Syringe Filters (Sterile)

Part Number	Description	Pkg
4902	MachV/0.2 µm Supor EKV membrane, 25 mm	50/pkg
4905	0.8/0.2 µm Supor membrane, 25 mm	50/pkg
4906	0.2 µm Ultipor membrane, 25 mm	50/pkg
4907	0.2 µm Fluorodyne II membrane, 25 mm	50/pkg
4908	0.2 µm Posidyne membrane, 25 mm	50/pkg

## Sterile Acrodisc Syringe Filters with Versapor Membrane



Ideal for clarification or prefiltration of virtually all aqueous samples.

- Variety of diameters and pore sizes to meet most aqueous sample filtration needs.
- Certified non-pyrogenic and biologically safe.

### Applications

- The economical choice for prefiltration of viscous or particulate-laden samples, serum filtration, and dissolution testing.

### Specifications

#### Materials of Construction

Filter Media: Versapor membrane (acrylic copolymer on a nonwoven support)  
Housing: Modified acrylic  
Effective Filtration Area  
25 mm: 2.8 cm<sup>2</sup>

#### Inlet/Outlet Connections

25 mm: Female luer lock inlet, male slip luer outlet

#### Typical Hold-up Volume

(with air purge)  
25 mm: > 50 µL

#### Maximum Operating Temperature

25 mm: 55 °C (131 °F)

#### Maximum Operating Pressure

25 mm: 5.2 bar (520 kPa, 75 psi)

#### Typical Water Flow Rate

mL/min at 3.1 bar (310 kPa, 45 psi)  
0.8 µm, 25 mm: 600  
1.2 µm, 25 mm: 660  
5 µm, 25 mm: 970

#### Endotoxin

< 0.25 EU/mL using Limulus Amoebocyte Lysate (LAL) test

#### Sterilization

Sterilized by gamma irradiation and individually packaged

### Ordering Information

Acrodisc Syringe Filters with Versapor Membrane (Sterile)

Part Number	Description	Pkg
4188	0.8 µm, 25 mm	50/pkg
4190	1.2 µm, 25 mm	50/pkg
4199	5 µm, 25 mm	50/pkg

## DMSO-Safe Sterile Acrodisc Syringe Filter with Nylon Membrane



Polypropylene housing and nylon membrane are compatible with solutions containing DMSO.

### Applications

- Ideal for sterilization of media used for cell cryopreservation.
- Offers broad chemical resistance to common solvents.

### Specifications

#### Materials of Construction

Filter Media: Nylon

Housing: Polypropylene

#### Effective Filtration Area

2.8 cm<sup>2</sup>

#### Inlet/Outlet Connections

Female luer lock inlet, male slip luer outlet

#### Typical Hold-up Volume

(with air purge)

≤ 50 µL

#### Maximum Operating Temperature

55 °C (131 °F)

#### Maximum Operating Pressure

6.2 bar (620 kPa, 90 psi)

#### Typical Water Flow Rate

60 mL/min at 2.1 bar (210 kPa, 30 psi)



#### Corporate Headquarters

Port Washington, NY, USA  
+1-800-717-7255 toll free (USA)  
+1-516-484-5400 phone

#### European Headquarters

Fribourg, Switzerland  
+41 (0)26 350 53 00 phone

#### Asia-Pacific Headquarters

Singapore  
+65 6389 6500 phone

#### Endotoxin

< 0.25 EU/mL using Limulus Amoebocyte Lysate (LAL) test

#### Sterilization

Sterilized by ethylene oxide and individually blister packed

### Ordering Information

#### DMSO-Safe Acrodisc Syringe Filter with Nylon Membrane (Sterile)

Part Number	Description	Pkg
4433	0.2 µm, 25 mm	50/pkg

### Sterile Acrodisc Syringe Filter Complementary Products

- Precise, high-performance centrifugal devices to process < 100 µL to 60 mL.
- Fast, sterile, AcroCap™ positive pressure device for volumes up to 3 liters.
- Efficient AcroPak™ 20 filters for volumes up to 5 liters.
- Innovative, sterile VacuCap® bottle-top filters for vacuum filtration of 100 mL to 5 liters.
- Capsules with Supor, Fluorodyne II, and Versapor membranes for filtration of larger volumes.

Pall Corporation has the most comprehensive family of scaleable filtration products that simplify validations and ease transitions from R&D to pilot to full-scale production. Look for these scaleable products with Supor, Supor EKV, Fluorodyne II, Posidyne, and Ultipor membranes:

- Membrane disc filters
- Acrodisc syringe filters with polypropylene housings
- AcroPak 20 filters for processing up to 5 liters
- AcroPak 200 capsules for processing 5 to 20 liters
- AcroPak 400 and 800 capsules for processing volumes over 10 liters
- Process-scale capsules and cartridges from Pall Corporation.

Visit us on the Web at [www.pall.com/laboratory](http://www.pall.com/laboratory)  
Contact us at [www.pall.com/contact](http://www.pall.com/contact)

Pall Corporation has offices and plants throughout the world. To locate the Pall office or distributor nearest you, visit [www.pall.com/contact](http://www.pall.com/contact).

The information provided in this literature was reviewed for accuracy at the time of publication. Product data may be subject to change without notice. For current information consult your local Pall distributor or contact Pall directly.

© Copyright 2022, Pall Corporation. Pall, , AcroCap, Acrodisc, AcroPak, Fluorodyne, Posidyne, Supor, Ultipor, VacuCap, and Versapor are trademarks of Pall Corporation.  
® Indicates a trademark registered in the USA.