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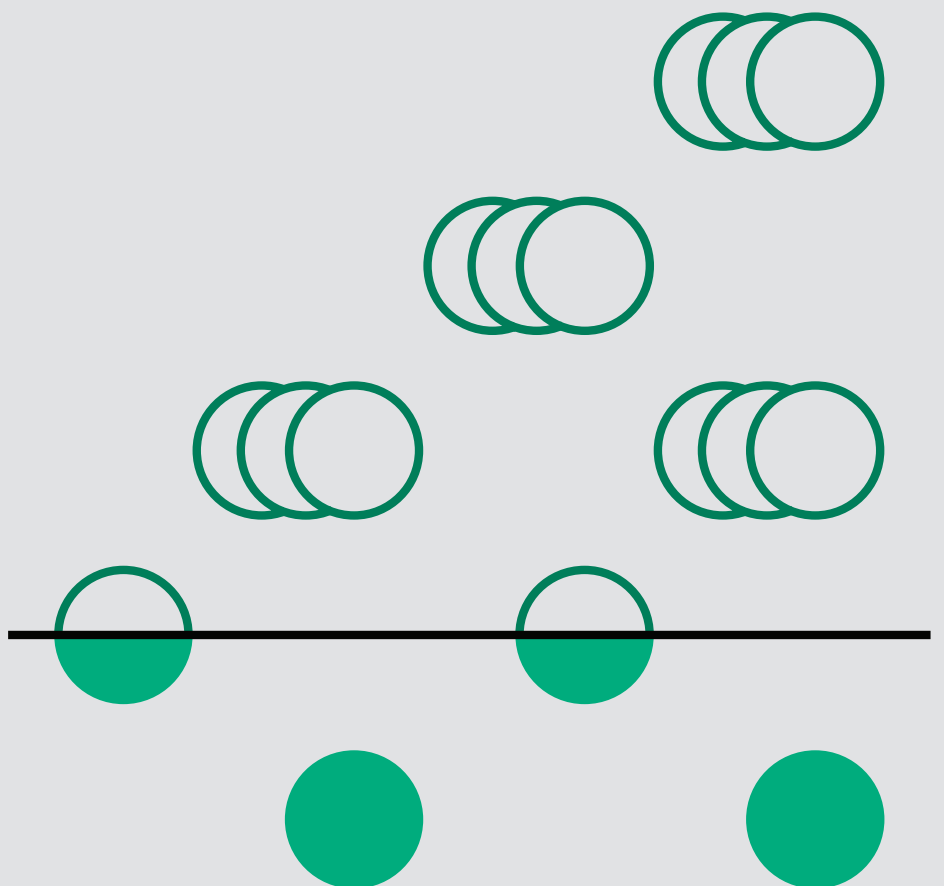
Bioprocess TFF(Tangential Flow Filtration)
solution cassette catalogue

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1	Ultrafiltration TFF	2
2	Microfiltration TFF	5
3	Single-Use TFF (SUTFF)	7
4	Single-Pass TFF (SPTFF): Inline-Concentrator	10
5	Single-Pass TFF (SPTFF): Modules	12
6	Single-Pass TFF (SPTFF): Inline-Diafiltration	16
7	TFF Holder Families	18

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Technical Information

	Omega™ (PES)	Delta (Cellulose)
Materials of Construction		
Membrane	Omega™ Polyethersulfone	Delta Regenerated Cellulose
Support	Polyolefin	
Screens	Polypropylene	
Encapsulant	Polyurethane with white pigment (TiO ₂)	
Permeate seals	Medical grade, platinum cured silicone	
Gaskets	Medical grade, platinum cured silicone	
Operating Limits		
Maximum pressure ⁽¹⁾	6 barg (87 psig) at 23 °C 4 barg (58 psig) at 55 °C	6 barg (87 psig) at 23°C 4 barg (58 psig) at 55°C
Maximum transmembrane pressure (TMP)	4 barg (58 psig) at 55 °C	4 barg (58 psig) at 55°C
Temperature range ⁽²⁾	-5 to 55 °C	4 to 55°C
pH range	2 to 14	2 to 13
Typical operating parameters		
Cross flow rate for processing	5 – 7 L/min/m ² (0.5 – 0.7 L/min/f ²)	5 to 7 L/min/m ² (0.5 to 0.7 L/min/ft ²)
Cross flow rate for cleaning	8 – 10 L/min/m ² (0.8 – 1.0 L/min/ft ²)	5 to 8 L/min/m ² (0.5 to 0.8 L/min/ft ²)
Integrity test		
Test pressure	2 barg (30 psig)	4.1 barg (60 psig) for 10, 30 kDa MWCO 2.1 barg (30 psig) for 100 kDa MWCO
Maximum air forward flow	< 1600 mL/min/m ² (< 150 mL/min/ft ²)	< 538 mL/min/m ² (< 50 mL/min/ft ²) for 10, 30 kDa MWCO ≤ 409 mL/min/m ² (≤ 38 mL/min/ft ²) for 100 kDa MWCO

⁽¹⁾ Pressure rating will be dependent on rating of the lowest system component.
⁽²⁾ Cassettes must not be allowed to freeze.

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Specification

	Omega™ (PES)	Delta (Cellulose)
Shelf Life		
Storage solution	0.3 N Sodium hydroxide	0.1 N Sodium hydroxide
Expected shelf life from	5 years	3 years
Storage condition	In the original packaging at 4-25°C and protected from direct light	

Biological Safety
The materials of construction in the T-Series cassettes have been tested and meet requirements for United States pharmacopeia (USP) biological reactivity test, in vivo <88> at 70°C
Documentation
Certificate of quality
Membrane Cassette Care and Use Procedures
Material safety data sheet (MSDS) for cassette preservative
Two platinum cured silicone gaskets

Ordering information

Membrane Type

Molecular Weight Cutoff (MWCO)

Nominal membrane area

T-Series cassettes

Part Number :

Example Part Number:

OS010T12

Omega™ membrane, 10 kDa molecular weight cutoff, 0.1 m² (1.1 ft²) Centramate™ screen channel cassette

Membrane Type		
Code	Membrane	Description
OS	Omega™	Modified Polyethersulfone
DC	Delta	Regenerated Cellulose

Molecular Weight Cutoff (MWCO)		
Code	NMWC	
001	1 kDa	
003	3 kDa	
005	5 kDa	
010 ⁽¹⁾	10 kDa	
030 ⁽¹⁾	30 kDa	
050	50 kDa	
070	70 kDa	
100 ⁽¹⁾	100 kDa	
300	300 kDa	

⁽¹⁾ Note: Delta membrane is only available in 10, 30, and 100 kDa MWCO.

Nominal membrane area		
Code	Cassette type	Nominal membrane area
T01	Centramate™	93 cm² (0.1 ft²)
T02	Centramate™	186 cm² (0.2 ft²)
T12	Centramate™	0.1 m² (1.1 ft²)
T06	Centrasette™	0.5 m² (5.4 ft²)
T26	Centrasette™	2.5 m² (27.0 ft²)

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Operating Limits	
Recommended maximum operating feed pressure	5 barg (75 psig)
Maximum operating temperature	50 °C (122 °F)

Cassette channel configuration	
Medium screen	Clarified solutions; low – medium viscosity Final concentration < 30% (300 mg/mL)
Suspended screen/open channel	Particles / cells / cell debris in solution / High viscosity

Storage agent	
Shipped wet in	15 – 20% glycerin, 0.05 – 0.1% sodium azide
Storage solution	0.3 N sodium hydroxide

The storage solution must be removed, and the cassette flushed well with water prior to use to prevent product contamination.

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PS

Example Part Number:
PSM10C12P2

Microfiltration cassettes with Supor™ membrane,
0.1 micron rating, medium screen,
0.02 m² (0.2 ft²) Centramate™ cassette

Membrane Type		
Code	Membrane	Description
PS	Supor™ PES	Microfiltration-optimized Polyethersulfone

Available micron ratings		
	Code	Pore size (µm)
	M10	0.1
	M20	0.2
	M45	0.45
	M65	0.65

Nominal membrane area		
Code	Cassette and screen type	Nominal membrane area
C12P1	Centramate™, Medium screen	0.01 m² (0.1 ft²)
C12P2	Centramate™, Medium screen	0.02 m² (0.2 ft²)
C12	Centramate™, Medium screen	0.1 m² (1.0 ft²)
F06	Centrasette™, Medium screen	0.5 m² (5.4 ft²)
F26	Centrasette™, Medium screen	2.5 m² (27.0 ft²)
C11P1	Centramate™, Suspended screen	0.01 m² (0.1 ft²)
C11P2	Centramate™, Suspended screen	0.02 m² (0.2 ft²)
C11	Centramate™, Suspended screen	0.1 m² (1.0 ft²)
F07	Centrasette™, Suspended screen	0.5 m² (5.4 ft²)
F27	Centrasette™, Suspended screen	2.5 m² (27.0 ft²)

Note: Not all varieties of PSM cassettes shown in the tables are available, it depends on size/micron rating.

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Technical Information

Omega™ (PES)	
Materials of Construction	
Membrane	Polyethersulfone
Support	Polyolefin
Screens	Polypropylene
Encapsulant	Polyurethane with white pigment (TiO₂)
Seals	Platinum cured silicone and thermoplastic elastomer
Manifold plates	Glass bead reinforced polypropylene with white pigment (TiO₂)
Port caps	Polypropylene for Luer connections, polyethylene for sanitary fittings
Covers	High impact polystyrene

Operating Limits	
Recommended maximum operating feed pressure	4 barg (58 psig)
Maximum transmembrane pressure (TMP)	4 barg (58 psig)
Temperature range during processing	4 to 40 °C
Temperature range during pre-use sanitization	18 to 25 °C
Recommended cross flow rate	5 – 7 L/min/m² (0.5 – 0.7 L/min/ft²)
Typical operating time	8 hours

Integrity test	
Test pressure	2 barg (30 psig)
Maximum air Forward Flow	≤ 1600 mL/min/m² (≤150 mL/min/ft²) at 20 °C

Single-use TFF modules must be used in an appropriate holder.
They are integrated in single-use TFF systems. The operating limits of the system should also be considered.

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CSUM

Example Part Number:
CSUM010T050

Single use TFF Module with omega membrane,
10 kDa nominal molecular weight cutoff,
0.5 m² (5.4 ft²) Centrasette™ screen channel cassette

Membrane Type		
Code	Membrane	Description
CSUM	Omega™	Single-use TFF Module with Omega™ membrane

Molecular Weight Cutoff (MWCO)		
Code	NMWC	
010	10 kDa	
030	30 kDa	
100	100 kDa	

Nominal membrane area		
Code	Cassette type	Nominal membrane area
T001	Centramate™	93 cm ² (0.1 ft ²)
T002	Centramate™	186 cm ² (0.2 ft ²)
T010	Centramate™	0.1 m ² (1.1 ft ²)
T050	Centrasette™	0.5 m ² (5.4 ft ²)
T250	Centrasette™	2.5 m ² (27.0 ft ²)

Ordering information - Holders

Holders for Single Use TFF Modules

For Centramate™ membrane type (0.1 m² membrane area)

Part Number :

CSUH040

Holder for Centramate™ type membrane
Orientation: Both (Vertical or horizontal)
Number of Single use TFF module that fit in holder: 1
Part numbers of compatible Single use TFF module:
CSUM__T001, CSUM__T002, CSUM__T010

For Centrasette™ membrane type (0.5 and 2.5 m² membrane area)

Part Number :

CSUH250SA

Holder for Centrasette™ type membrane
Orientation: Horizontal
Number of Single use TFF module that fit in holder: 1
Part numbers of compatible Single use TFF module:
CSUM__T050, CSUM__T250

Cadence™ Inline Concentrator

Ultrafiltration, Ready for use, SPTFF

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Technical Information

Materials of Construction	
Cassette	Omega™ (PES) or Delta (Regenerated Cellulose) membrane
Screen	Polypropylene
Encapsulant	Polyurethane with white pigment (TiO ₂)
Permeate seals	Medical grade silicone
Gaskets	Medical grade, platinum-cured silicone
Manifold plates	Ultra high molecular weight polyethylene (UHMWPE)
Tubing and connectors	Tubing: pharmaceutical grade platinum-cured silicone T01/T02 - female Luer connection and cap: polypropylene (PP) T12 - MPC quick disconnect connector and plug: polysulfone (PS) T06 - MPX quick disconnect connector and plug: polysulfone (PS)
Operating Limits	
Maximum operating pressure	4.1 barg (60 psig)
Maximum transmembrane pressure (TMP)	4.1 barg (60 psig)
Operating temperature range	20 – 25 °C for extended use (freezing will damage module) Up to 40 °C for 4 hours cleaning
Operating pH range	2 to 13
Integrity test	
Delta membrane module	< 538 mL/min/m ² (<50 mL/min/ft ²) at 4.1 barg (60 psig)
Omega™ membrane module	<1600 mL/min/m ² (<150 mL/min/ft ²) at 2.1 barg (30 psig)

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Example Part Number:

ILD010T120407

Cadence™ Inline concentrator cassette with Delta regenerated cellulose membrane, 10 kDa molecular weight cutoff, 0.7 m² (7.5 ft²) membrane area

Membrane Type		
Code	Membrane	Description
ILOS	Omega™	Inline concentrator cassette™ with Omega™ PES membrane
ILD	Delta	Inline concentrator cassette™ with Delta regenerated cellulose membrane

Molecular Weight Cutoff (MWCO)	
Code	NMWC
010	10 kDa
030	30 kDa

Nominal membrane area	
Code	Membrane area
T010	0.065 m ² (0.7 ft ²)
T020	0.13 m ² (1.4 ft ²)
T120	0.7 m ² (7.5 ft ²)
T060	3.5 m ² (37.7 ft ²)

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
Technical Information

Delta (Cellulose)	
Materials of Construction	
Membrane	Regenerated Cellulose
Gaskets	USP Class VI Pharmaceutical grade, platinum cured silicone
Manifold plates	USP Class VI Ultra-high-molecular-weight (UHMW) polyethylene
Connections	1⁄8 inch hose barb
Operating Limits	
Feed pressure during processing	2.1 to 4.1 barg (30 to 60 psig)
Retentate pressure during processing	Greater than 0.3 barg (5 psig)
Maximum operating pressure	6 barg (87 psig) at 23 °C 4 barg (58 psig) at 55 °C
Operating temperature range	4 to 40 °C (122 °F)
Operating pH range	2 to 13
Integrity test	
Test pressure	4.1 barg (60 psig)
Maximum air Forward Flow	< 500 mL/min/m² (<50 mL/min/ft²) at 20 °C
Storage agent	
Storage solution	0.1 N sodium hydroxide for less than three months 5% ethanol and 1% sodium diacetate for more than three months

Format guide

Module Area (m²)	Cadence™ Single Pass TFF Configurations			
	4-in-series	7-in-series	8-in-series	9-in-series
T01 (93 cm²)	0.065	0.12	0.14	0.17
T02 (186 cm²)	0.13	0.24	0.28	0.33
T12 (0.1 m²)	0.7	1.3	1.5	1.8
T06 (0.5 m²)	3.5	6.5	7.5	9.0

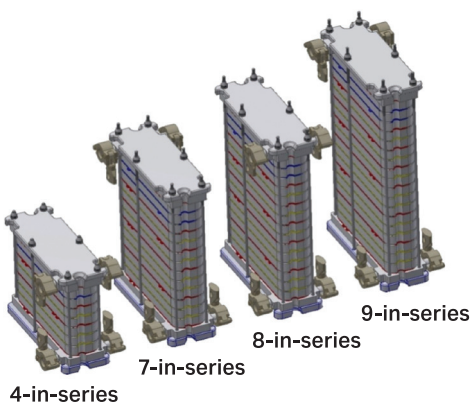Modules shown utilizing T12 Centramate™ (0.1 m²) building blocks

Ordering information

Membrane Type	Molecular Weight Cutoff (MWCO)	Nominal membrane area
---------------	--------------------------------	-----------------------

Cadence™ SPTFF Modules

Part Number :

Example Part Number:

CD010T120713

Cadence™ Single pass TFF Module with Delta membrane, 10 kDa molecular weight cutoff, 1.3 m² (14.0 ft²) membrane area

Membrane Type		
Code	Membrane	Description
CD	Delta	Cadence™ Single-pass TFF Module with Delta membrane

Molecular Weight Cutoff (MWCO)		
	Code	NMWC
	010	10 kDa
	030	30 kDa

Nominal membrane area and configuration		
Code	Configuration	Membrane area
T010407	4-in-series	0.065 m ² (0.7 ft ²)
T010713	7-in-series	0.12 m ² (1.3 ft ²)
T010815	8-in-series	0.14 m ² (1.5 ft ²)
T010918	9-in-series	0.17 m ² (1.8 ft ²)
T020407	4-in-series	0.13 m ² (1.4 ft ²)
T020713	7-in-series	0.24 m ² (2.6 ft ²)
T020815	8-in-series	0.28 m ² (3.0 ft ²)
T020918	9-in-series	0.33 m ² (3.6 ft ²)
T120407	4-in-series	0.13 m ² (1.4 ft ²)
T020713	7-in-series	0.24 m ² (2.6 ft ²)
T020815	8-in-series	0.28 m ² (3.0 ft ²)
T020918	9-in-series	0.33 m ² (3.6 ft ²)
T120407	4-in-series	0.7 m ² (7.5 ft ²)
T120713	7-in-series	1.3 m ² (14.0 ft ²)
T120815	8-in-series	1.5 m ² (16.1 ft ²)
T120918	9-in-series	1.8 m ² (19.4 ft ²)

Ordering information - Holders

Holders for Cadence™ SPTFF Module

Part Number :

CSPH

Example Part Number:

CSPHS9000

Cadence™ Single-Pass Holder for Centrasette™ T06 format
Compatible with module configuration : 7-in-series, 8-in-series, 9-in-series


Holder Type	
Code	Description
CSPH	Cadence™ Single-Pass Holder

Cassette Format and Maximum area		
Code	Cassette Format	Maximum Membrane Area
M0330	Centramate™ (T01, T02)	0.33 m ² (3.6 ft ²)
M1800	Centramate™ (T12)	1.8 m ² (19.4 ft ²)
S3500	Centrasette™ (T06)	3.5 m ² (37.7 ft ²)
S9000	Centrasette™ (T06)	9.0 m ² (96.9 ft ²)

Application

Single-Pass TFF Modular Kit

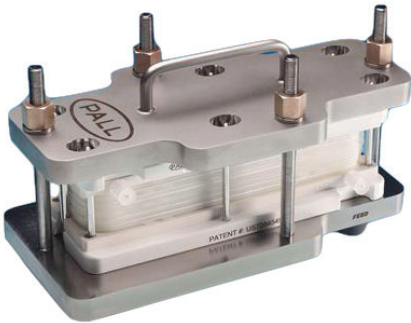- 93 cm² or 186 cm² nominal area cassettes can be used
- Process Development scale Only
- Flexibility to create any number of stages and staging formats up to 9-in-series
- 1 Cadence™ modular kit, 1 Holder, TFF cassette needed

Ordering information

Cadence™ Modular Kit Part Number :

SPTFFMKPD9

Kit Contents: 1 x Feed manifold
 1 x Retentate manifold
 25 x Gaskets


Holder Part Number (ordered separately) :

CSPHM0330

Cadence™ SPTFF holder

FS001K10

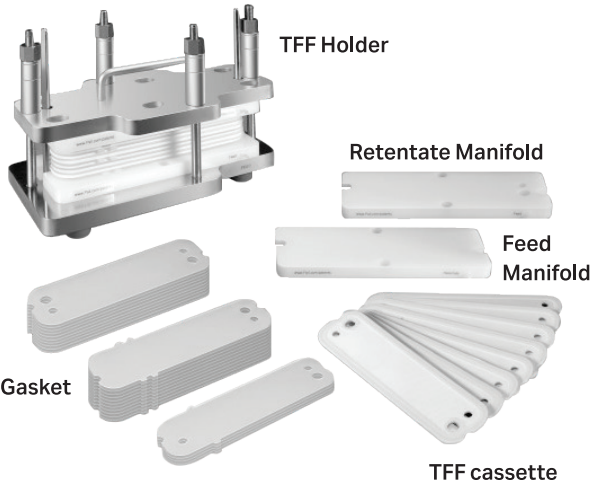Centramate™ standard TFF holder

Number of cassettes each holder type can accommodate:

	Cadence™ SPTFF Holder		Centramate™ Holder	
	Minimum	Maximum	Minimum	Maximum
T01	4	30	1	21
T02	3	21	1	15

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Technical Information

Materials of Construction	
Cassette	Omega™ (PES) or Delta (Regenerated Cellulose) membrane
Screen	Polypropylene
Encapsulant	Polyurethane with white pigment (TiO ₂)
Permeate seals	Medical grade silicone
Gaskets	Medical grade, platinum-cured silicone
Manifold plates	Ultra high molecular weight polyethylene (UHMWPE)
Connections	T01/T02 – female Luer connection and cap: Polypropylene T12 – female MPC quick disconnect connector and plug: Polysulfone
Flowpaths	Polypropylene, polysulfone, and platinum-cured silicone
Pump Head Tubing	PharmaPure (platinum-cured silicone) or Sta-Pure (expanded polytetrafluoroethylene (ePTFE) and platinum-cured silicone) Flowpath connections on tubing sets are open to the feed container/vessel
Operating Limits	
Maximum operating pressure	4.1 barg (60 psig)
Operating temperature range	15 – 25 °C for extended use (freezing may damage the module) Up to 40 °C for 4 hours cleaning
Operating pH range	2 to 13
Integrity test	
Delta membrane module	< 538 mL/min/m ² (<50 mL/min/ft ²) at 4.1 barg (60 psig)
Omega™ membrane module	<1600 mL/min/m ² (<150 mL/min/ft ²) at 2.1 barg (30 psig)

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Membrane Type		
Code	Membrane	Description
DFOS	Omega™	Inline diafiltration cassette with Omega™ PES membrane
DFDC	Delta	Inline diafiltration cassette with Delta regenerated cellulose membrane

Molecular Weight Cutoff (MWCO)	
Code	NMWC
030	30 kDa

Nominal membrane area	
Code	Membrane area
T010	0.11 m ² (1.2 ft ²)
T020	0.22 m ² (2.4 ft ²)
T120	1.2 m ² (12.9 ft ²)

TFF Holder Families

Centramate™, Centrasette™, Centrastak™

Lab and Development Scale (Centramate™)

	Centramate™ LV	Centramate™ SS	Centramate™ PE
Maximum Operating Temperature	50 °C	80 °C	80 °C
Maximum Operating Pressure	5 bar	6 bar	6 bar
Surface Finish	Ra <0.8µm - Passivated	Ra <0.8µm - Passivated	Ra <0.8µm - Passivated
Nominal Membrane Area	0.01 – 0.04 m²	0.01 – 0.37 m²	0.01 – 0.46 m²
Dimensions (mm)	230(L) x 120(W) x 66(H)	230(L) x 135(W) x 160(H)	240(L) x 95(W) x 170(H)
Working Volume (starting)	25 – 4,000 mL	0.5 – 80 L	0.5 – 100 L
Working Volume (Final)	15 mL	150 mL	150 mL
Weight	4 kg	6.5 kg	7 kg

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cytiva.co.kr

Cytiva and the Drop logo are trademarks of Life Sciences IP Holdings Corporation or an affiliate. Cadence, Centramate, Centrasette, Contrastak and Supor are trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva. Any third-party trade marks are the property of their respective owners.

© 2024 Cytiva.

For local office contact information, visit cytiva.com/contact

CY47806-10Dec24-BR