


Monitor pH/C-900


Short Instructions

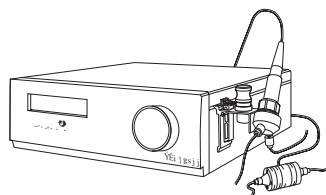

Step	Action
1	<p>Switch on the instrument by means of the mains switch on the rear panel.</p> <div>Calibrating</div>
2	<p>The main operating menu (RUN-menu) is shown.</p> <div> Cond Temp pH 25.4% 22.9C 11.5 </div>
3	<p>Calibrate the pH electrode before use and/or daily by using 2 buffers with known pH values.</p> <div>Calibrate pH</div>
4	<p>The conductivity flow cell does not normally need to be calibrated.</p>

Step	Action
5	Set the conductivity scaling for 0 and 100%. <div>Set Cond Scale 100%</div>
6	Always store the pH electrode in 1:1 mixture of pH 4 buffer and 1 M KNO ₃ when not in use.

Menu overview Monitor pH/C-900

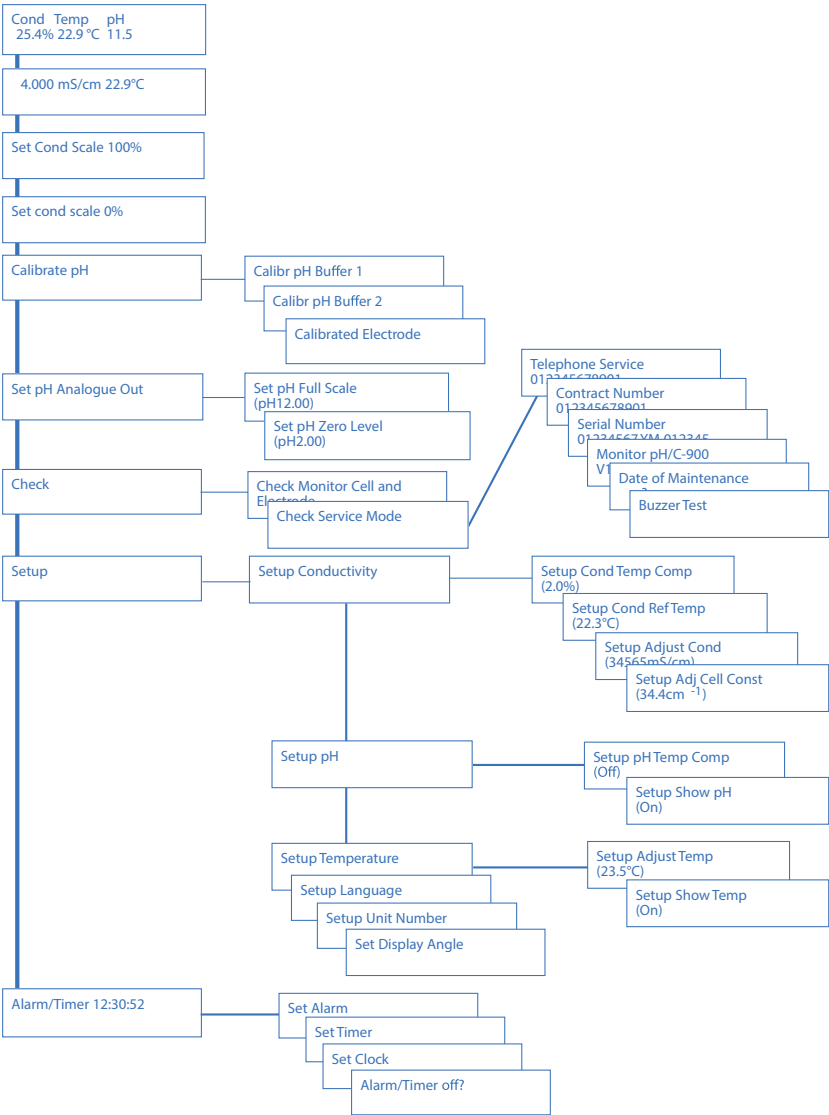



cytiva.com

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

All other third-party trademarks are the property of their respective owners.

© 2020–2021 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

18112010 AF V:5 06/2021