


Mounting kit for Pump P-950/P-960


Instruction

CAUTION! Before installing Pump P-950 or Pump P-960 in an ÄKTAdesign™ system, contact your local GE Healthcare Bio-Sciences service organisation to make sure that the system is equipped with the correct electrical components.


Installation

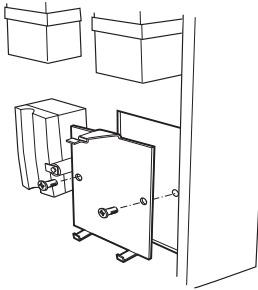


1. Attach the mounting plate to the lower right-hand corner at the inside of the system door, using the two flat-headed screws.

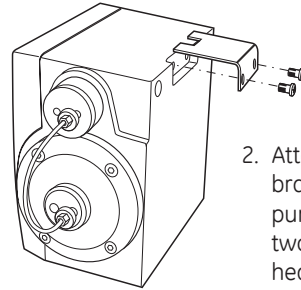

2. Attach the plate bracket to the pump using the two fillister-head screws.

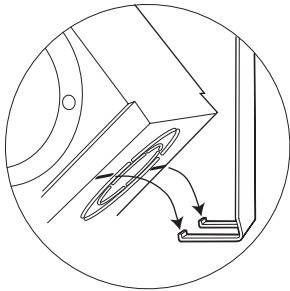


3. Fit the two hooks on the mounting plate in the holes underneath the pump.

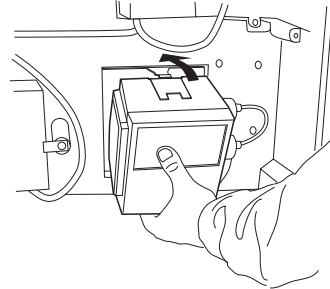

4. Push the pump towards the mounting plate as the figure shows. A click sound will indicate that the pump is locked in position.

Removing the pump

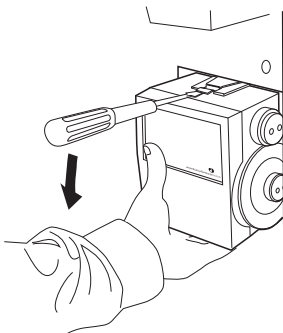

1. Release the locking plate with a screw driver while holding the pump.
2. Remove the pump.



WARNING! PINCH HAZARD

Do not grip the pump with the fingers at the rear of the pump.



www.gehealthcare.com
GE Healthcare Bio-Sciences AB
Björkgatan 30
751 84 Uppsala
Sweden

ÄKTAdesign is a trademark of GE Healthcare Ltd, a General Electric company. GE, imagination at work and GE monogram are trademarks of General Electric Company.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. GE Healthcare reserves the right, subject to any regulatory and contractual approval, if required, to make changes in specifications and features shown herein, or discontinue the product described at any time without notice or obligation. Contact your local GE Healthcare representative for the most current information.

© 2005 General Electric Company - All rights reserved.

GE Healthcare Bio-Sciences AB, a General Electric company.

GE Healthcare
Munzinger Strasse 5, D-79111 Freiburg, Germany

GE Healthcare
Amersham Place, Little Chalfont, Buckinghamshire, HP7 9NA, UK

GE Healthcare
800 Centennial Avenue, P.O. Box 1327, Piscataway, NJ 08855-1327, USA

GE Healthcare
Sanken Bldg. 3-25-1, Hyakunincho, Shinjuku-ku, Tokyo 169-0073, Japan



imagination at work