

# Tubing Connector

(Code No. 19-2150-01)

## Connection of capillary tubing

1. Cut the tubing cleanly at a 45° angle.
2. Remove the screw nut and slip it over the tubing.
3. Remove the sealing plug.
4. Push the tubing through the sealing plug until it protrudes about 1 cm.
5. Seat the tubing in the nipple. Slide the sealing plug along the tubing and into the nipple.
6. Finger-tighten the connector.



[www.gehealthcare.com](http://www.gehealthcare.com)

GE Healthcare Bio-Sciences AB  
Björkgatan 30  
751 84 Uppsala  
Sweden

GE Healthcare Europe GmbH  
Munzinger Strasse 5  
D-79111 Freiburg  
Germany

GE Healthcare UK Ltd  
Amersham Place  
Little Chalfont  
Buckinghamshire, HP7 9NA  
UK

GE Healthcare Bio-Sciences Corp  
800 Centennial Avenue  
P.O. Box 1327  
Piscataway, NJ 08855-1327  
USA

GE Healthcare Bio-Sciences KK  
Sanken Bldg.  
3-25-1, Hyakunincho  
Shinjuku-ku, Tokyo 169-0073  
Japan

GE, imagination at work and GE monogram are trademarks of General Electric Company.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. GE Healthcare reserves the right, subject to any regulatory and contractual approval, if required, to make changes in specifications and features shown herein, or discontinue the product described at any time without notice or obligation. Contact your local GE Healthcare representative for the most current information.

© 2006 General Electric Company - All rights reserved.

GE Healthcare AB, a General Electric Company.



imagination at work

56-4745-00 AE 05/2006