

Monitor UVis-920

Appendix to User Manual, BioProcess system control



1 Appendix

1.1 Monitor UVis-920, BioProcess system control

Monitor UVis-920 is a UV and visible light absorption monitor for use in liquid chromatography. The instrument is designed as a stand-alone instrument but can as option be equipped with a process card and an extended rear panel for use in a BioProcess™ system.

The process card offers possibilities to control UVis-920 with digital signals or an external display box, and receive UV absorbance data on a 4-20 mA output.

The User Manual 11-0031-12 is valid also for the BioProcess system control version of the UVis-920. This appendix gives additional information only.

1.2 Connecting electrical cables

The sockets for electrical signals are located on the rear panel.


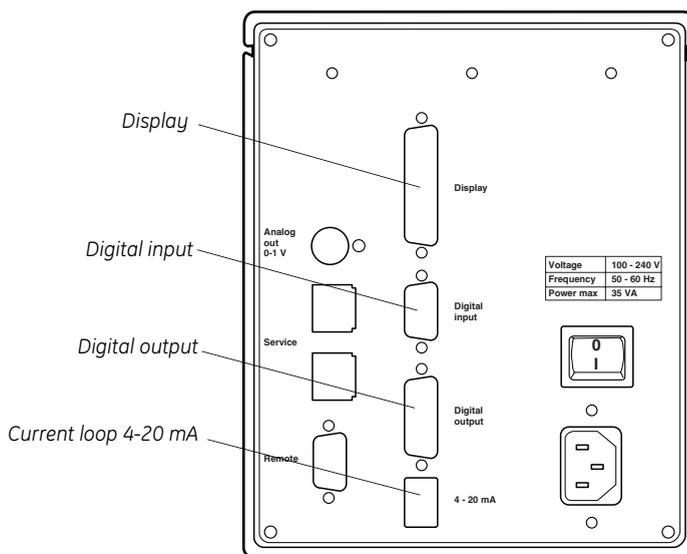


Fig 1-1. UVis-920 rear panel for use in a BioProcess system

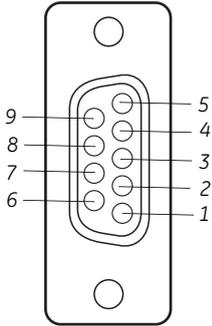

1.2.1 Connecting digital input signals


Connect digital input signals to the 9-pole D-Sub male **Digital input** connector. All digital inputs are galvanic isolated from UVis-920 main electronic. All inputs uses TTL-levels by default. With jumper W1 on the process card the alter level can be changed to 10 volt.

Pin	Signal	Function
1	Set range bit 0	Set AU-range for 4-20 mA output
2	Set range bit 1	
3	Set range bit 2	
4	Set range bit 3	
5	Auto zero	0=autozero
6	Run/End	0=run, 1=end
7	Keyboard lock	0=display locked 1=display works normally
8	Voltage 5 or 12 V (max 100 mA)	Output voltage for external pull-ups
9	Ground 0 V	

Table 1-1.

1.2.2 Connecting digital output signals

All digital outputs are placed in the 15-pole D-Sub female **Digital output** connector and functions as a relay that closes to ground. Hence external pull-ups are needed. All digital outputs are galvanic isolated from UVis-920 main electronic.

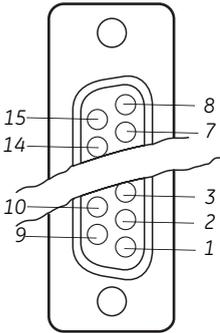

Pin	Signal	Function
1	Get range bit 0	Shows AU-range for 4-20 mA
2	Get range bit 1	
3	Get range bit 2	
4	Get range bit 3	
5	Filter code bit 0	Shows code for inserted filter
6	Filter code bit 1	
7	Filter code bit 2	
8	Filter code bit 3	
9	Error code on display	0=error, 1=no error
10	Run end	0=running, 1=end state
11	Out of range	0=AU-value is out of range 1=AU-value is in range
12	Ground 0 V	
13		
14	24 V max 100 mA	Output voltage for external pull-ups
15		

Table 1-2.

1.2.3 Binary range codes for digital inputs and outputs

Binary value range code				Range
Bit 3	Bit 2	Bit 1	Bit 0	
0	0	0	0	Not used (when this value is set on digital input, range can be changed from display)
0	0	0	1	(-1 6) AU
0	0	1	0	(-0.5 3) AU
0	0	1	1	(-0.25 1.5) AU
0	1	0	0	(-0.125 0.75) AU
0	1	0	1	(-1 1) AU
0	1	1	0	(0 1) AU
1	1	1	1	Not used (when this value is set on digital input, range can be changed from display)

Table 1-3. Binary range codes for digital inputs and outputs


1.2.4 Binary filter codes for digital outputs

Binary filter code				Filter wavelength
Bit 3	Bit 2	Bit 1	Bit 0	
0	0	0	0	Filter not fitted
1	0	0	1	546 nm
1	0	1	0	405 nm
1	0	1	1	280 nm
1	1	0	0	260 nm
1	1	0	1	215 nm
1	1	1	0	Custom filter
1	1	1	1	No filter inserted

Table 1-4. Binary filter codes for digital outputs

1.2.5 Connecting equipment to 4-20 mA output

Connect any equipment to the 3-pole female connector **Current loop 4-20 mA**.

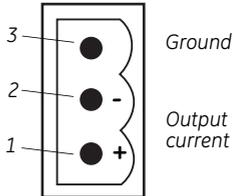

Pin	Signal	Function
1	+	Output current
2	-	Output current
3	GRD	Ground

Table 1-5.


1.3 Error messages

Error codes 201-206 are internal errors that the BioProcess interface shows on its display.

Message	Description/Action
Error 201 Not connected	BioProcess interface has no connection to UVis-920. Call GE Healthcare.
Error 202 Internal error	Internal error. Call GE Healthcare.
Error 203 Communication	Communication protocol between BioProcess and UVis-920 differ. Call GE Healthcare.
Error 204 Communication	BioProcess interface has no connection to UVis-920. Call GE Healthcare.
Error 205 Communication	BioProcess could not connect to UVis-920 at start-up. Call GE Healthcare.
Error 206 Serial no not valid	Flow Cell serial number entered from the BioProcess display is not valid

Table 1-6.

1.4 Menu overview, BioProcess system control

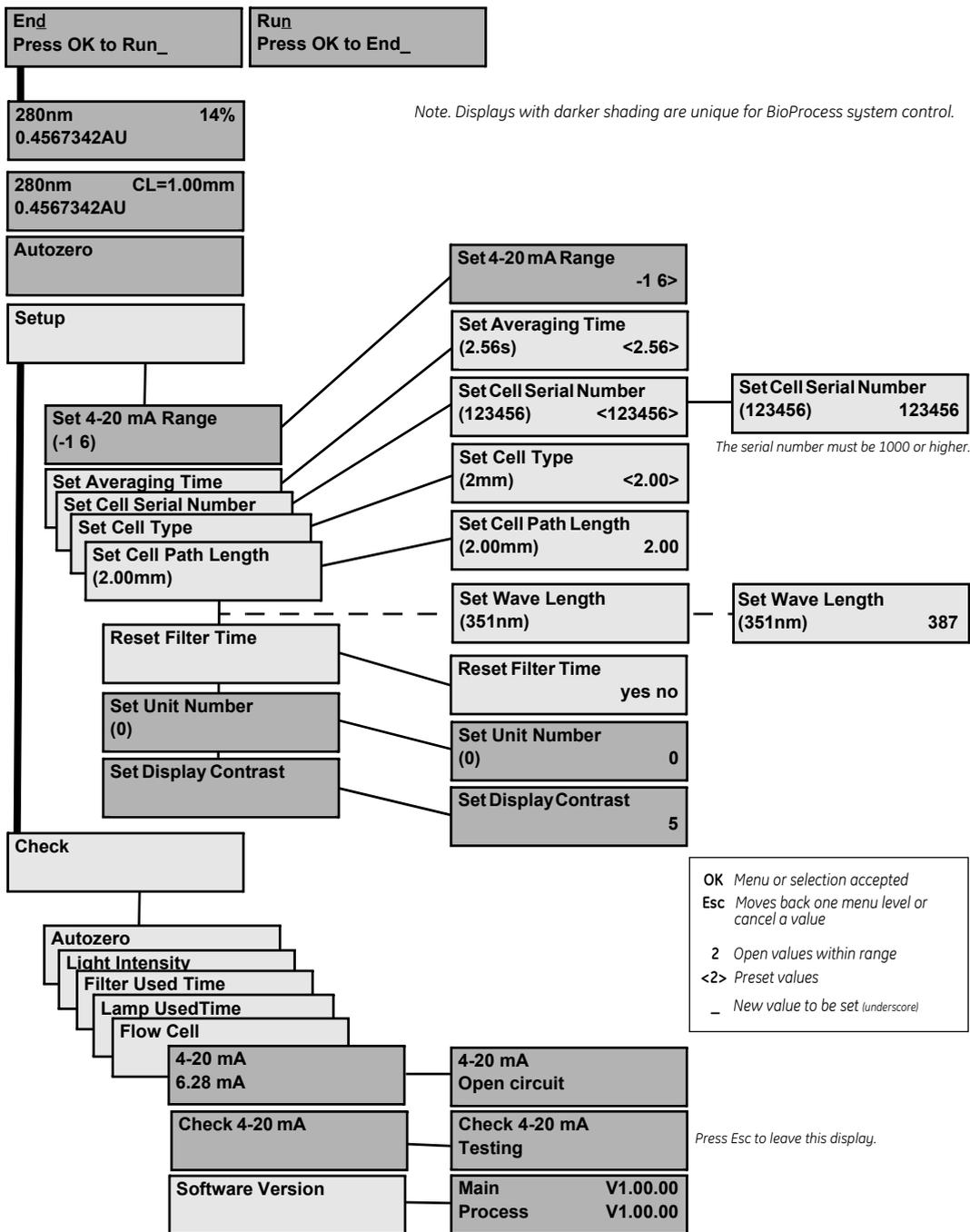

1.5 Menus

Most of the windows (messages) that the BioProcess interface shows on its display are identical with the windows presented on the stand alone UVis-920 display, except the one marked with darker shading in the menu overview. These are unique for the BioProcess system control interface. See 1.4 Menu overview, BioProcess system control.

1.5.1 Main menu

To start and stop the monitor (compare with the run/end key on the monitor):

End
Press OK to Run_

1 Start measuring (run mode), press **OK**.

Run
Press OK to End_

2 Stop measuring (end mode), press **OK**.

280nm **14%**
0.4567342AU

The display shows the filter wavelength in nm and the absorbance value in AU. The percentage of the set 4-20 mA range that the absorbance value corresponds to is also indicated.

Example: If 4-20 mA is set to (0 1) and the absorbance value is 0.5 AU, the display shows 50%.

280nm **CL=1.00mm**
0.4567342AU

Indicates the filter wavelength in nm, cell path length in mm and the current AU value.

Autozero

The autozero display/function sets the detected absorbance to zero when **OK** is pressed.

1.5.2 Setup menu

Set 4-20 mA range

This menu allows you to select range for the 4-20 mA output.

- 1 Select menu **Setup** and press **OK**.
- 2 Select menu **Set 4-20 mA Range**, press **OK**.
- 3 Set the range in AU, press **OK**.

Set 4-20 mA Range
(-1 6)

Set 4-20 mA Range
-1 6>

Value	Range
-1 6	-1 to +6.0 AU
-0.5 3	-0.5 to + 3.0 AU
-0.25 1.5	-0.25 to +1.5 AU
-0.125 0.75	-0.125 to +0.75 AU
-1 1	-1.0 to +1.0 AU
0 1	0 to +1.0 AU

Set unit number

The unit number for UVis-920 can be selected by the user.

- 1 Select menu **Setup** and press **OK**.
- 2 Select menu **Set Unit Number**, press **OK**.
- 3 Set unit number, press **OK**. Values: 0-9.

Set Unit Number
(0)

Set Unit Number
(0) <9>

Set display contrast

The display contrast can be selected by the user.

- 1 Select menu **Setup**, press **OK**.
- 2 Select menu **Set Display Contrast**, press **OK**.
- 3 Set contrast, press **OK**. Values: 0-9.

Set Display Contrast

Set Display Contrast
5

1.5.3 Check menu**Indication of 4-20 mA output**

The current value of the 4-20 mA output can be indicated.

- 1 Select menu **Check**, press **OK**.
- 2 Select sub menu **4-20 mA**, press **OK**. The display shows the current value in mA.

4-20 mA
6.28 mA

4-20 mA
Open circuit

If the 4-20 mA output is not used (not connected) the display shows **Open circuit**.

Check 4-20 mA output


The function of the 4-20 mA output can be tested.

- 1 Select menu **Check**, press **OK**.
- 2 Select sub menu **Check 4-20 mA**, press **OK**.
- 3 Start the test, press **OK**. The test will ramp the output up to 20 mA and then decrease the signal in 10% steps back to 4 mA. The test is run continuously.

Check 4-20 mA

Check 4-20 mA
Testing

Stop the test by pressing **Esc**.

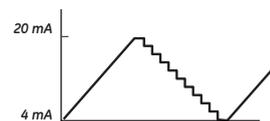

Software version

- 1 Select menu **Check**, press **OK**.
- 2 Select menu **Software Version**, press **OK**. The display shows the software version of the main system and the process system.

Main	V1.00.00
Process	V1.00.00

www.chromatography.amershambiosciences.com

GE Healthcare
Amersham Biosciences AB
Björkgatan 30
751 84 Uppsala
Sweden

Drop Design and BioProcess are trademarks of GE Healthcare Ltd.
GE and GE monogram are trademarks of General Electric Company.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. GE Healthcare reserves the right, subject to any regulatory and contractual approval, if required, to make changes in specifications and features shown herein, or discontinue the product described at any time without notice or obligation. Contact your local GE Healthcare representative for the most current information.

© 2005 General Electric Company – All rights reserved.

Amersham Biosciences AB, a General Electric company going to market as GE Healthcare.

GE Healthcare Amersham Biosciences AB
Björkgatan 30, 751 84 Uppsala, Sweden

GE Healthcare Amersham Biosciences Europe GmbH
Munzinger Strasse 9, D-79111 Freiburg, Germany

GE Healthcare Amersham Biosciences UK Ltd
Amersham Place, Little Chalfont, Buckinghamshire, HP7 9NA, UK

GE Healthcare Amersham Biosciences Corp
800 Centennial Avenue, P.O. Box 1327, Piscataway, NJ 08855-1327, USA

GE Healthcare Amersham Biosciences KK
Sanken Bldg. 3-25-1, Hyakunincho, Shinjuku-ku, Tokyo 169-0073, Japan

Asia Pacific Tel: +852 2811 8693 Fax: +852 2811 5251 • Australasia Tel: + 61 2 9899 0999 Fax: +61 2 9899 7511 • Austria Tel: 01/57606-1619 Fax: 01/57606-1627 • Belgium Tel: 0800 73 888
Fax: 03 272 1637 • Canada Tel: 800 463 5800 Fax: 800 567 1008 • Central, East, & South East Europe Tel: +43 1 982 3826 Fax: +43 1 985 8327 • Denmark Tel: 45 16 2400 Fax: 45 16 2424 • Finland &
Baltics Tel: +358-(0)9-512 39 40 Fax: +358 (0)9 512 39 439 • France Tel: 01 69 35 67 00 Fax: 01 69 41 96 77 • Germany Tel: 0761/4903-490 Fax: 0761/4903-405 • Italy Tel: 02 27322 1 Fax: 02 27302 212
• Japan Tel: +81 3 5331 9336 Fax: +81 3 5331 9370 • Latin America Tel: +55 11 3933 7300 Fax: +55 11 3933 7304 • Middle East & Africa Tel: +30 210 9600 687 Fax: +30 210 9600 693 • Netherlands
Tel: 0165 580 410 Fax: 0165 580 401 • Norway Tel: 815 65 555 Fax: 815 65 666 • Portugal Tel: 21 417 7035 Fax: 21 417 3184 • Russia & other C.I.S. & N.I.S Tel: +7 (095) 232 0250, 956 1137 Fax: +7 (095)
230 6377 • South East Asia Tel: 60 3 8024 2080 Fax: 60 3 8024 2090 • Spain Tel: 93 594 49 30 Fax: 93 594 49 55 • Sweden Tel: 018 612 1900 Fax: 018 612 1910 • Switzerland Tel: 0848 8028 12 Fax:
0848 8028 13 • UK Tel: 0800 616928 Fax: 0800 616927 • USA Tel: 800 526 3593 Fax: 877 295 8102



imagination at work