

# ÄKTExpress large column holder

## Installation instructions

Step	Action
1	Use the enclosed hex key to loosen the two screws on the LEFT column holder of ÄKTExpress system approximately half a turn.
2	Remove the column holder from the system by pulling it to the right.
3	Mount the large column holder by pushing it into the correct position on the LEFT side of the system.
4	Fasten the large column holder by tightening the two screws with the hex key.

**WARNING**

The large column holder must **ONLY** be mounted on the **LEFT** side of the system. Do **NOT** mount it on the right side of the system as the large column will hinder the function of the fraction collector.



**CORRECT.** The large column holder correctly mounted.



**WRONG.** Do NOT mount the large column holder on the right side of the system!

Page intentionally left blank



## cytiva.com

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

ÅKTA™ is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

All other third-party trademarks are the property of their respective owners.

© 2020 – 2021 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit [cytiva.com/contact](https://cytiva.com/contact)

28400974 AB V:4 05/2021