turvaohjeet

Kasetti

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(Suomi)

Miten kasettipussi avataan turvallisesti

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(English)

How to Open the Cassette Bag Safely

Go to www.gehealthcare.com for complete handling and storage instructions in other languages

The cassette and cassette bag contain an aqueous solution containing up to 0.2 N NaOH and 21% glycerin. When opening the cassette bag, follow the standard safety procedures for handling aqueous NaOH, including the use of safety glasses, safety gloves, and protective lab coat.

To avoid unexpected spilling, hold the cassette package upright over a sink and trim the top of the cassette bag with scissors. Drain and dispose of the excess storage solution in accordance with environmental regulations.

To avoid inadvertent contact with the storage solution after the cassette

is removed from the bag, rinse the outside surface of the cassette with high purity water.

If you allow an ultrafiltration cassette to dry out, the membrane will be damaged. Therefore, do not store the cassette without rewetting it with an approved storage solution. Read the handling and storage instructions in the guide entitled: *GE Healthcare Cassettes: Quick Reference Guide*.

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