


Pump P-905

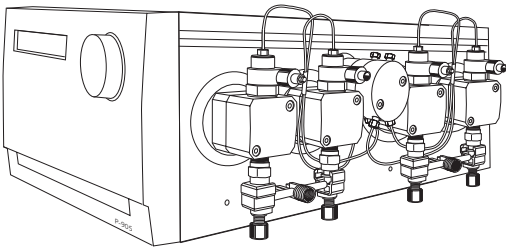


- 1 **Switch on the instrument** by means of the mains switch on the rear panel.
- 2 Press **OK** to start synchronisation.
- 3 The **main operating menu** is shown.
- 4 Set the **flow rate**.
- 5 Start the pump by setting it in **RUN**-mode.
- 6 The **main operating menu** is shown.
- 7 Start a gradient by setting the **Gradient Length** and **Gradient Target**.
- 8 Stop the pump by setting it in **END**-mode.

Selftest Please wait...	
Pump synchronisation Press OK to start	
End	0.00ml/min 0.00 MPa 0.0%B
Set Flow Rate (0.000ml/min) 0.5	
End	0.500ml/min Run
Run	2.000ml/min 2.00 MPa 50.0%B
Set Gradient Length (0.00min) 8.0	
Set Gradient Target (0.0%B) 20.0	
Run	0.500 ml/min End Hold Pause



Menu overview Pump P-905

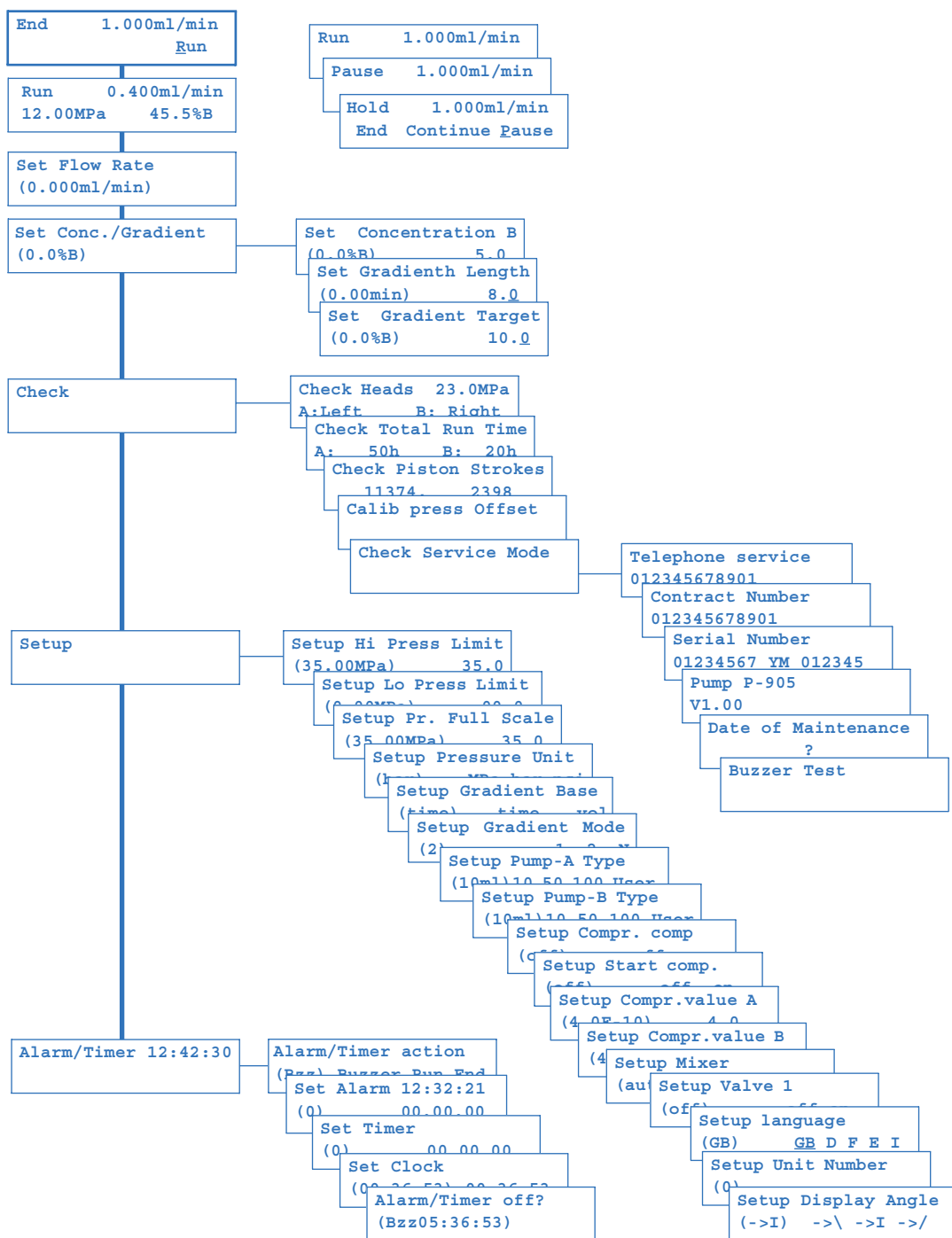

GE imagination at work

GE Healthcare Bio-Sciences AB, a General Electric Company.
GE Healthcare Bio-Sciences AB
Björkgatan 30
751 84 Uppsala, Sweden
© 2006 - 2009 General Electric Company—All rights reserved.