


Monitor UV-900

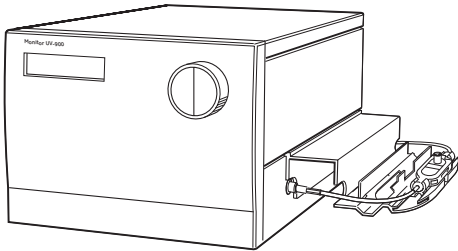

Action


- 1 **Switch on the module** with the mains switch on the rear panel.
- 2 **Switch on the UV-lamp** by selecting **Run**.
- 3 **Select wavelength** for one, two or three wavelengths.
- 4 **Select Autozero** and press **OK**.
- 5 **The main operating menu** (RUN-menu) is shown.

Display

Calibrating Please wait...
End Press OK to Run
Set Wavelength (215, 254, 280 nm)
Set Wavelength λ 1 (215 nm)
Set Wavelength λ 2 (254 nm)
Set Wavelength λ 3 (280 nm)
Autozero
λ 1 (215) 1.12345 AU λ 2 (254) 0.02345 AU



Menu overview Monitor UV-900

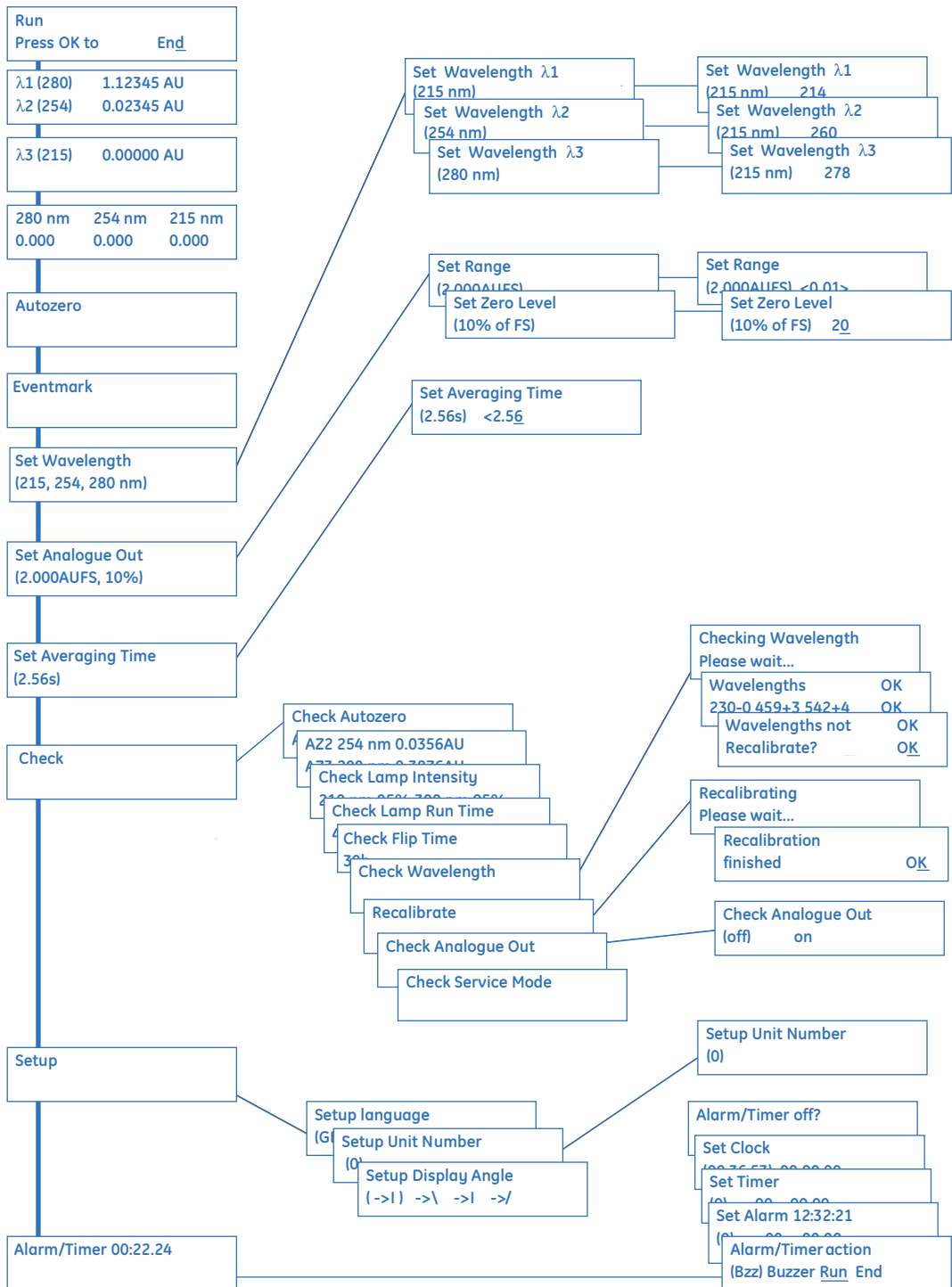

GE imagination at work

GE Healthcare Bio-Sciences AB, a General Electric Company.
GE Healthcare Bio-Sciences AB
Björkgatan 30
751 84 Uppsala, Sweden
© 2006 - 2009 General Electric Company—All rights reserved.