


Monitor pH/C-900

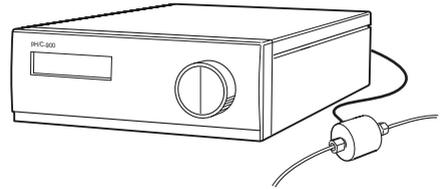

Action

- 1 **Switch on the instrument** by means of the mains switch on the rear panel.
- 2 The **main operating menu** (RUN-menu) is shown.
- 3 The conductivity flow cell does not normally need to be calibrated.
- 4 Set the **conductivity scaling** for 0 and 100%.

Display


Calibrating
Please wait...

Cond	Temp
25.4%	22.9 °C

Set Cond Scale 100%
(50.00 mS/cm)



Menu overview Monitor pH/C-900

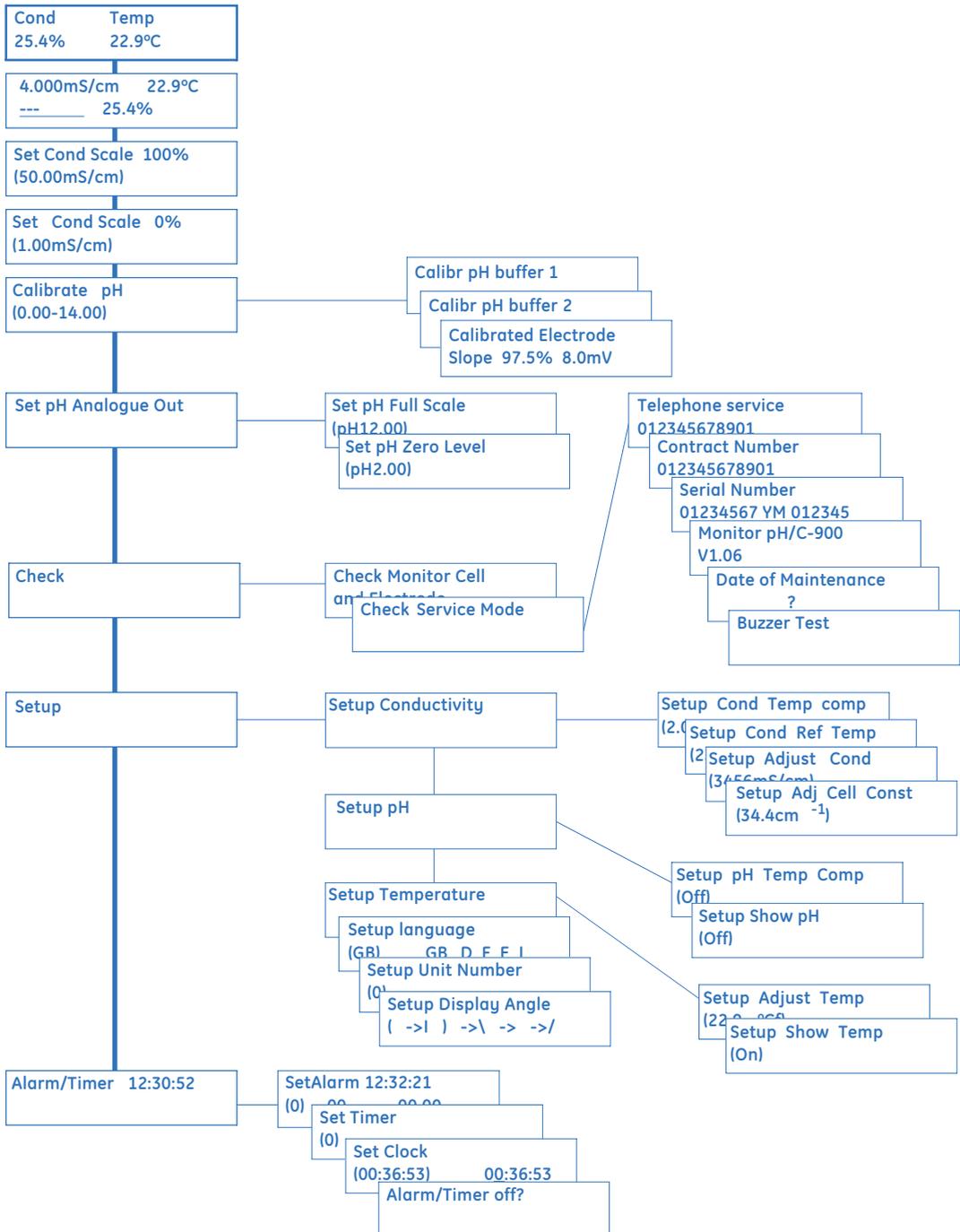

GE imagination at work

GE Healthcare Bio-Sciences AB, a General Electric Company.
 GE Healthcare Bio-Sciences AB
 Björkgatan 30
 751 84 Uppsala, Sweden
 © 2006 - 2009 General Electric Company—All rights reserved.