


ReadyClamp


Instructions

ReadyClamp is a single use clamp designed to fit the ReadyMate™ Disposable Aseptic Connector (DAC) and other 1½" sanitary fittings that conform to ISO Standard 2852 for Sanitary Couplings. ReadyClamp is a light-weight clamp for traceable, tamper proof connections.


| Step | Action |
|------|--|
| 1 | Remove the clamp from the packaging and open it. If needed, record the serial number printed on the cable tie. Notice the location marker on the cable tie shown as "XXX". |

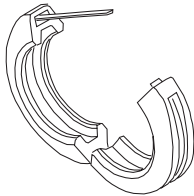


| | |
|---|--|
| 2 | Fit the clamp around DAC connection. Align the frosted area and the arrow of the DAC toward the opening of the clamp. Use hand pressure to wrap the clamp around the DAC and properly seat the DAC in the clamp channel. Hold the clamp in position. |
|---|--|

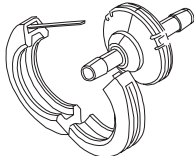

| Step | Action |
|------|---|
| 3 | While holding the DAC in position, pull back the free end of the black cable tie and thread it into the black latching head of the tie located in the opposing clamp. Pull on the cable tie end to secure the clamp around the DAC. |

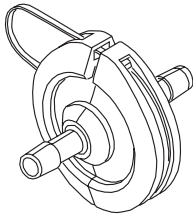

| | |
|---|--|
| 4 | Use a cable tie tensioning tool to tighten the cable tie until at least one "X" on the strap becomes visible. Once this point is achieved, apply tension and twist cable tie tensioning tool to cut the strap. The connection is now complete. |
|---|--|

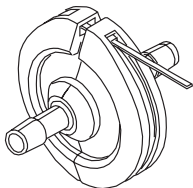

Note: Apply maximum 50 pounds of cable tension.

Technical specifications

| | |
|-----------------------|--|
| Outside clamp profile | 50 mm (1 ½"), ISO standard 2852 |
| Materials | Clamp: Polycarbonate Cable tie: Polyamide |

Associated documentation

| | |
|---|----------|
| ReadyMate Disposable Aseptic Connector (DAC) Instructions | 28938570 |
|---|----------|

Ordering information

| | Quantity | Code no. |
|----------------|----------|----------|
| ReadyClamp 1½" | 25 | 28936690 |



cytiva.com

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

ReadyToProcess and ReadyMate are trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

All other third-party trademarks are the property of their respective owners.

© 2020–2021 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

28940620 AB V:5 02/2021