

# Modification note for Injection valve INV-917

(Code No. 18-1147-23)

## **Please check this after removal of the distributing plate:**

If the valve's switching plate is 9 mm thick, the valve housing has to be modified for correct spring pressure for the switching plate of Valve Kit 18-1147-56 (10 mm thick).

Please contact GE Healthcare service (Labcrew™) to perform the following:

1. Remove the valve housing (see valve service manual 18-1116-41).
2. Note how the shaft is orientated vis-à-vis the ports, then push out the shaft and **remove two of the three 0.5 mm washers.**
3. Refit.

Labcrew is a trademark of GE Healthcare companies. GE, imagination at work and GE monogram are trademarks of General Electric Company.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. Contact your local GE Healthcare representative for the most current information.

© 2006 General Electric Company - All rights reserved.

GE Healthcare Bio-Sciences AB, a General Electric Company.

GE Healthcare Bio-Sciences AB, Björkgatan 30, 751 84 Uppsala, Sweden



imagination at work