

Prefilter

The prefilter (Product Code 19508401) is used to filter buffers and solvents from the pump to the column at pressure up to 10 MPa (100 bar, 1420 psi).



Table of contents

Unpacking	3
Connecting the prefilter	3
Replacing the filter	3
Spare parts	3

1. Unpacking

Please check the contents of the package against this list.

Designation	No. supplied
Prefilter (see Fig 1.)	1
Preflanged capillary tubing, ETFE	1
Instructions	1

2. Connecting the prefilter

Connect the prefilter between the pump and the injection valve in the direction shown in Figure 1.

3. Replacing the filter

1. Unscrew the two sections of the prefilter.
2. Remove the O-ring and filter with a laboratory spatula.
3. Place a new filter into the prefilter's female part and press it down with the spatula.
4. Press the O-ring into place, sealing the edges all the way around the filter.
5. Screw the two parts of the prefilter back together and reconnect the prefilter.

4. Spare parts

Designation	Code No.	No. per pack
Filters + O-rings (prefilter)	19-5082-01	5+2

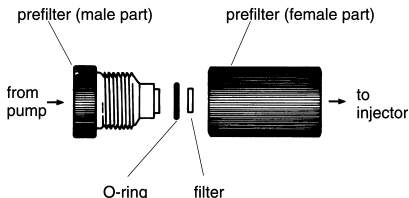


Fig 1. Positioning and component parts of the prefilter.

For local office contact
information, visit,
www.gelifesciences.com/contact

GE Healthcare Bio-Sciences AB
Björkgatan 30
751 84 Uppsala
Sweden

www.gelifesciences.com

GE Healthcare Europe GmbH
Munzinger Strasse 5
D-79111 Freiburg
Germany

GE Healthcare UK Ltd
Amersham Place
Little Chalfont
Buckinghamshire, HP7 9NA
UK

GE Healthcare Bio-Sciences Corp
100 Results Way, Marlborough
MA 01752
USA

GE Healthcare Dharmacon, Inc.
2650 Crescent Dr., Lafayette
CO 80026
USA

HyClone Laboratories, Inc.
925 W 1800 S, Logan
UT 84321
USA

GE Healthcare Japan Corporation
Sanken Bldg. 3-25-1, Hyakunincho
Shinjuku-ku, Tokyo 169-0073
Japan

GE and GE monogram are trademarks of General Electric Company.

© 1993-2016 General Electric Company
Previously published Oct. 1993

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.

