

M3 Screw

Before connecting SuperFrac™ to the LC Controller 2252, the screws in the communication cable, Code No. 19-6005-02, must be replaced.

Take one end of the communication cable, Code No. 19-6005-02, and replace the mounted screws with the M3 screws supplied. Connect that end to LC Controller 2252.

The M3 screws cannot be used for connection of the other GE Healthcare instruments.



For local office contact information, visit
www.gelifesciences.com/contact

GE Healthcare Bio-Sciences AB
Björkgatan 30
751 84 Uppsala
Sweden

www.gelifesciences.com

GE Healthcare Europe GmbH
Munzinger Strasse 5,
D-79111 Freiburg,
Germany

GE Healthcare UK Ltd
Amersham Place
Little Chalfont
Buckinghamshire, HP7 9NA
UK

GE Healthcare Bio-Sciences Corp
800 Centennial Avenue
P.O. Box 1327 Piscataway,
NJ 08855-1327
USA

GE Healthcare Japan corporation
Sanken Bldg.
3-25-1, Hyakunincho
Shinjuku-ku, Tokyo 169-0073
Japan

GE, imagination at work and GE monogram are trademarks of General Electric Company.
SuperFrac is a trademark of GE Healthcare companies.

© 1993-2010 General Electric Company – All rights reserved.
Previously published Apr. 1993.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.



imagination at work

56-1061-05 AH 09/2010