

# Attach Rail extension and Column holders

## Instructions

### Scope

Instructions on how to attach the following Rail extension and column holders to the instrument.


**Note:** *The described column holders are designed for the ÄKTA™ avant and ÄKTA pure instruments and may not be possible to attach to other ÄKTA instruments. Also, column holders used for other ÄKTA instruments may not be possible to attach to the ÄKTA avant and ÄKTA pure instruments.*

Holder	Illustration
Rail extension (29011352)	
Column holder rod (28956270)	

Holder	Illustration
Column holder (28956282)	
Column clamp o.d. 10-21 mm (28956319)	

## Attach Rail extension and Column holder rod

Press down the button on the rod (1) and attach the rod to the rails of the instrument (2).

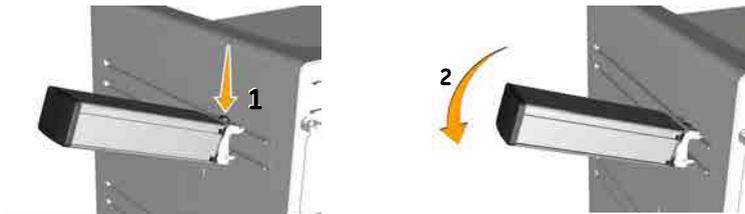


# Attach Column holder

Step	Action
------	--------

- |   |  |
|---|--|
| 1 | Attach the lower hooks of the column holder snap-in to the lower rail. |
|---|--|



- |   |  |
|---|--|
| 2 | Press the tab on the holder and attach the upper hooks of the column holder snap-in to the upper rail. |
|---|--|

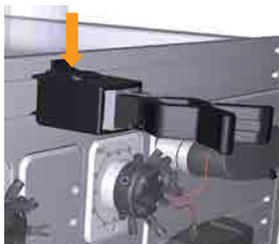

# Attach Column clamp

Step	Action
------	--------

- |   |   |
|---|---|
| 1 | Press the inner end tabs of the clamp together. |
|---|---|



Step	Action
------	--------

---

- 2 Insert the clamp into a rail of the instrument.



- 3 Turn the clamp 90 degrees and release.



Page intentionally left blank



## cytiva.com/akta

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

AKTA is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

All other third-party trademarks are the property of their respective owners.

© 2020 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit [cytiva.com/contact](https://cytiva.com/contact)

28968641 AD V:4 11/2020