

ÄKTA avant and ÄKTA pure

Attach sample cassettes to the instrument

Instructions

Scope

Instructions on how to attach cassettes to the ÄKTA™ avant and ÄKTA pure instruments. All cassettes designed for the fraction collector can be attached and used as, for example, sample containers.

Attach cassettes to the ÄKTA avant instrument



To install a cassette follow the instructions in the table below.

Step	Action
------	--------

- | | |
|---|---|
| 1 | Locate the four hooks below the instrument display. |
|---|---|



Step	Action
2	<p>Cassettes can be attached with either the short or long side facing the instrument.</p> <p>Hook on a cassette, placing one of the large side holes over two of the hooks.</p>

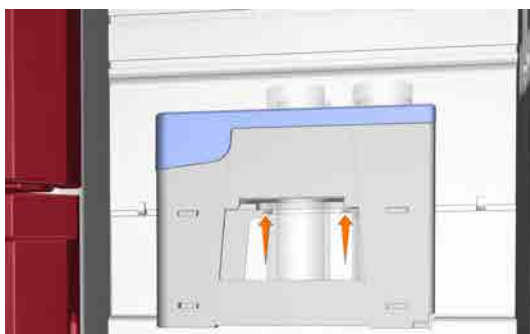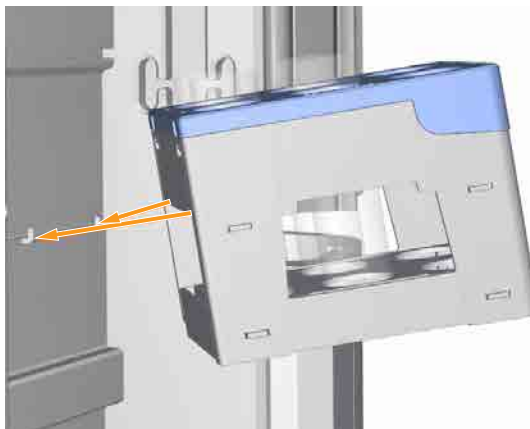


Attach cassettes to the ÄKTA pure instrument


To install a cassette follow the instructions in the table below. The Multi-purpose holder (29011349) must be used to attach a cassette to the ÄKTA pure instrument, see *Attach Multi-purpose holder and Loop holder (29021464)* for more information.

Step	Action
------	--------

- | | |
|---|---|
| 1 | Snap the upper hooks on the Multi-purpose holder into the upper holder rail on the left side of the instrument. |
|---|---|


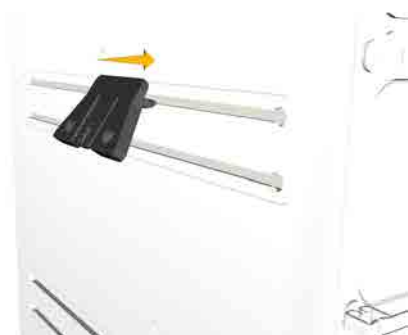


- | | |
|---|---|
| 2 | Press the tab on the holder and snap the lower hooks of the Multi-purpose holder into the lower rail. |
|---|---|

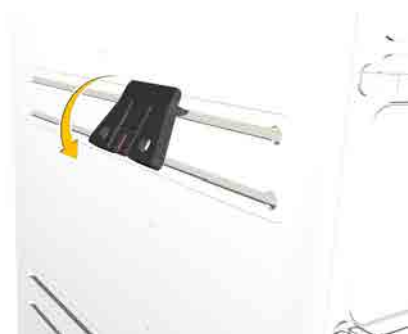

Step	Action
------	--------

- | | |
|---|---|
| 3 | <p>Cassettes can be attached with either the short or long side facing the instrument.</p> <p>Hook on a cassette, placing one of the large side holes over the top of the Multi-purpose holder.</p> |
|---|---|

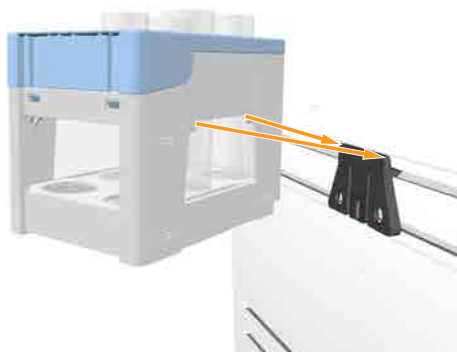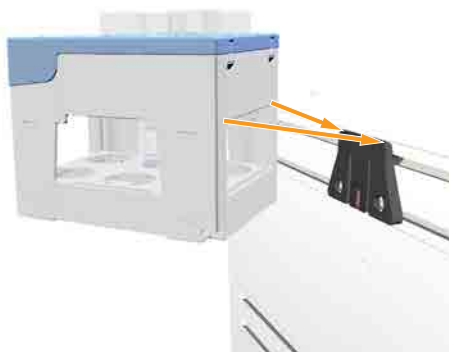

Page intentionally left blank



cytiva.com/akta

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

AKTA is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

All other third-party trademarks are the property of their respective owners.

© 2020–2021 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

28969192 AC V:5 01/2021