

Attach Multi-purpose holder and Loop holder

Instructions

Scope

Instructions on how to attach Multi-purpose holder (29011349) and Loop holder (29011350).


| Holder | Illustration | Used with... |
|------------------------------------|---|------------------------------|
| Multi-purpose holder (29011349) |  | ÄKTA™ avant and ÄKTA pure |
| Loop holder (29011350) |  | ÄKTA pure |

Description

Multi-purpose holder

The Multi-purpose holder can be used to attach accessories, eg., a Loop holder, tubing holders or a cassette. Attach the holder to a holder rail.

The illustration below shows the Multi-purpose holder.

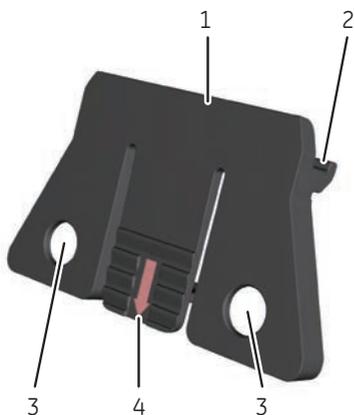

| Part | Function |
|------|--------------------------------------|
| 1 | Attachment point for accessories |
| 2 | Snap-in to holder rails |
| 3 | Attachment points for tubing holders |
| 4 | Tab |

Loop holder

The Loop holder can be used to attach up to five 10 ml capillary loops. Use two Multi-purpose holders to attach the holder to a holder rail.

The illustration below shows the Loop holder.



| Part | Function |
|------|--|
| 1 | Upper attachment to multi-purpose holder |
| 2 | Lower attachment to multi-purpose holder |

Installation

Attach Multi-purpose holder

Attach the holder to the rails of the instrument as described in the instructions below.


| Step | Action |
|------|--------|
|------|--------|

- | | |
|---|---|
| 1 | Snap the upper hooks on the Multi-purpose holder into the upper rail. |
|---|---|



| Step | Action |
|------|--------|
|------|--------|

- | | |
|---|---|
| 2 | Press the tab on the holder and snap the lower hooks of the Multi-purpose holder into the lower rail. |
|---|---|

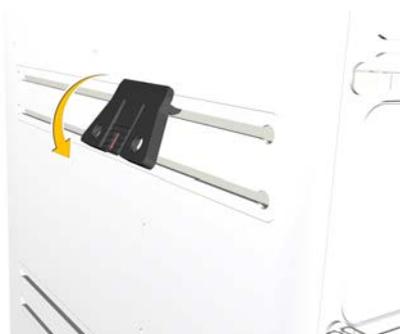

Result:

The Multi-purpose holder is attached.



Attach Loop holder

Attach the Loop holder to an ÄKTA pure instrument as described in the instructions below.

| Step | Action |
|------|--------|
|------|--------|

- | | |
|---|--|
| 1 | Attach two Multi-purpose holders to the rails, one on top of the other, within the area indicated below. |
|---|--|



| Step | Action |
|------|--------|
|------|--------|

- | | |
|---|---|
| 2 | Attach the hooks on the Loop holder to the two Multi-purpose holders, as illustrated below. |
|---|---|



Result:

The Loop holder is attached.





cytiva.com/akta

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

AKTA is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

All other third-party trademarks are the property of their respective owners.

© 2020–2021 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

29021464 AC V:5 01/2021