

Depletion of albumin from serum samples using ÄKTA start

Intellectual Property Notice: The Biopharma business of GE Healthcare was acquired by Danaher on 31 March 2020 and now operates under the Cytiva[™] brand. Certain collateral materials (such as application notes, scientific posters, and white papers) were created prior to the Danaher acquisition and contain various GE owned trademarks and font designs. In order to maintain the familiarity of those materials for long-serving customers and to preserve the integrity of those scientific documents, those GE owned trademarks and font designs remain in place, it being specifically acknowledged by Danaher and the Cytiva business that GE owns such GE trademarks and font designs.

cytiva.com

GE and the GE Monogram are trademarks of General Electric Company.

Other trademarks listed as being owned by General Electric Company contained in materials that pre-date the Danaher acquisition and relate to products within Cytiva's portfolio are now trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate. All other third-party trademarks are the property of their respective owners. © 2020 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

CY12781-20Apr20-AN

Application note 29-0642-95 AA


Depletion of albumin from serum samples using ÄKTA™ start


Albumin comprises 55% to 65% of the total protein in serum. The abundance of this protein makes it difficult to study other serum proteins, hence removal of albumin in a single step will speed up the process and aid in easy identification and characterization of low-abundance serum proteins, such as disease-specific biomarkers. Yang used UNICORN[™] start 1.0 control software to develop a method for quick and easy removal of albumin from human serum using ÄKTA start chromatography system, Frac30 fraction collector and a C 10/10 column packed with Blue Sepharose[™] 6 Fast Flow (FF) medium (resin).

Introduction

Serum is comprised of thousands of proteins and the composition varies with different physiological or pathological conditions. Some serum proteins may be potential biomarkers that are used for monitoring either disease progression or treatment. Effective depletion of albumin is a key step in the isolation and analysis of these biomarkers, which are typically present in low amounts.

Yang is a research scholar in a research laboratory focusing on identifying and characterizing disease-specific biomarkers in human serum. Yang felt too much of his time was being spent on manual purification tasks such as continuously monitoring the run at each step, collecting fraction manually and identifying protein fractions in the absence of an on-line UV detection facility. To save time and to make the process simple and effective, Yang switched to ÄKTA start with a Frac30 fraction collector. He used UNICORN start software from a neighboring lab to develop a method to remove albumin from human serum. Blue Sepharose is a chromatography medium commonly used to selectively remove albumin from serum samples or culture media. Blue Sepharose 6 FF is Cibacron™ Blue 3G coupled to Sepharose 6 FF. Albumin binds to Cibacron Blue 3G, a synthetic polycyclic dye that acts as an aromatic anionic ligand binding the albumin via electrostatic and/or hydrophobic interactions. Albumin depletion in turn enriches the other serum proteins, which are usually present only in very low concentrations.

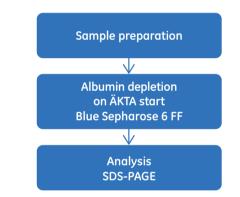

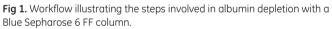

Methods

Yang followed the workflow illustrated in Figure 1 to perform albumin depletion using ÄKTA start. The depletion method was developed using the UNICORN start Method Editor module. The optimized run parameters Yang used are described in Table 1.



Table 1. UNICORN start method overview

| Method flow | Method settings |
|---------------------------|--|
| Method settings | User defined column, Column volume: 3 ml Pressure: 0.1 MPa ¹ Flow rate: 3 ml/min |
| Prime and equilibration | 3 CV ² |
| Sample application | Apply sample using loop, Sample volume: 1 ml Flow rate: 0.5 ml/min Fixed volume fractionation, Fraction volume: 1 ml |
| Wash out unbound | 3 CV ² Flow rate: 1 ml/min Fixed volume fractionation, Fraction volume: 1 ml |
| Elution and fractionation | Isocratic elution: % B concentration 100%, Elution volume: 5 CV ² Fractionation; Peak fractionation, level based (start/end: 10 mAU) |
| Prime and equilibration | 3 CV ² |

² 0.5 MPd = 5 bdf (45.5 ps
² CV = Column volumes

Yang exported the method from UNICORN start to a USB memory stick and imported the method into ÄKTA start instrument placed in a cold cabinet. Yang manually packed a C 10/10 column (i.d. 10 mm) with Blue Sepharose 6 FF medium (3 ml). The column was mounted onto ÄKTA start instrument and equilibrated with binding buffer (20 mM sodium phosphate buffer, pH 7.0). A sample loop was used to load 1 ml of sample (human serum) onto the column. The sample was loaded at a flow rate of 0.5 ml/min to ensure efficient binding of albumin to the media. Flowthrough was collected as 1 ml fixed volume fractions using Frac30 fraction collector. To remove unbound proteins, the column was washed with 3 column volumes (CV) of binding buffer at 1 ml/min flow rate.

Prepare ÄKTA start system by priming with desired buffers

Bound albumin was eluted from the column using 5 CV of 100% elution buffer (20 mM sodium phosphate buffer, 2 M NaCl, pH 7.0) in a single step. The protein peak was fractionated using UV level-based fractionation and 1.5 ml fractions were collected using Frac30 fraction collector. The run was monitored in real time using the UV trace from the instrument display. Using the bitmap (.bmp) result file, Yang identified the protein fractions collected during the sample application, wash unbound and elution steps. The pooled fractions were analyzed by 12.5% SDS polyacrylamide gel electrophoresis (SDS-PAGE).

Results

In this study Yang depleted albumin from serum samples in a single depletion step, using a C 10/10 column manually packed with Blue Sepharose 6 FF media. Yang took advantage of the additional options available in UNICORN start to develop a method that included multiphase fractionation. Fixed volume fractionation was used to collect flowthrough and peak fractionation (UV level-based) was used to collect bound albumin. The compact size of ÄKTA start, and the ability to run the instrument in a standalone mode (i.e. without being connected to UNICORN start software) with the Frac30 fraction collector allowed Yang to perform the depletion run in a cold cabinet. The Frac30 fraction collector allowed him to collect fractions at multiple phases without manual intervention.

Customize your method using UNICORN start Method Editor

To quickly view the results and to identify the peak fractions for pooling Yang printed the bitmap (.bmp) image file that had been generated during the purification run and saved onto a USB memory stick (Fig 2). He also imported the result file from the USB memory stick into UNICORN start (Fig 3A) and used Evaluation module functions such as peak integration for further analysis of the results and PDF report generation for project archival.

The SDS-PAGE profile of serum fractions showed > 95% removal of albumin from the human serum sample (Fig 3B). The flowthrough was free from albumin and suitable for further characterization.



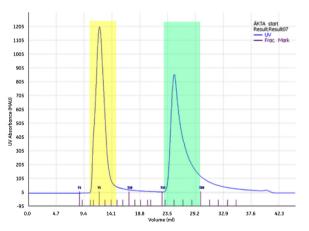


Fig 2. Chromatographic profile (.bmp) of albumin depletion from human serum sample on ÄKTA start. The highlighted areas (yellow: Flowthrough; green: Eluate) represent the pooled fractions.

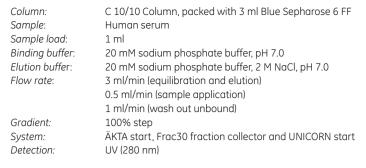

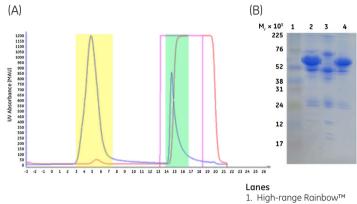

Ordering information

| Product | Quantity | Code number |
|--|------------|-------------|
| ÄKTA start | 1 | 29-0220-94 |
| UNICORN start 1.0 software | 1 | 29-0187-51 |
| Frac30 fraction collector | 1 | 29-0230-51 |
| Blue Sepharose 6 Fast Flow | 50 ml | 17-0948-01 |
| C 10/10 column | 1 | 19-5001-01 |
| High-range Rainbow molecular weight markers | 1 × 250 µl | RPN756E |

Related literature

| Product | Code number |
|---|-------------|
| Purification of N-terminal histidine-tagged protein using ÄKTA start, Application note | 29-0642-77 |
| Purification of GST-tagged protein using ÄKTA start, Application note | 29-0642-98 |
| Purification of antibodies using ÄKTA start and HiTrap Protein G HP column, Application note | 29-0643-02 |

Fig3. (A) Chromatogram showing albumin depletion from human serum sample on ÄKTA start (data imported from USB memory stick to UNICORN start software). The area highlighted in yellow represents the flowthrough fractions, and the green-highlighted area is the albumin-containing eluate). The red lines represents the conductivity curve and the pink line shows the gradient concentration. (B) SDS-12.5% PAGE profile of the sample following albumin depletion.

molecular weight markers

2. Load 3 Flowt

Flowthrough
Eluate

Summary

Yang used the intuitive UNICORN start software to quickly develop a method for effective depletion of albumin from human serum using ÄKTA start, Frac30 fraction collector, and a manually-packed column with Blue Sepharose 6 FF media. By using an automated process, Yang no longer needed to spend so much time monitoring the run. The bitmap result file generated by ÄKTA start enabled quick and easy identification of relevant fractions, and automatic fraction collection with Frac30 saved Yang time and ensured that the relevant fractions could be easily pooled.

Depletion of albumin helped Yang enrich the serum sample, facilitating biomarker isolation and characterization. Yang is now able to purify multiple batches to obtain low-abundance serum proteins for further characterization.

For local office contact information, visit **www.gelifesciences.com/contact**

www.gelifesciences.com/AKTAstart

GE Healthcare Bio-Sciences AB Björkgatan 30 751 84 Uppsala Sweden



imagination at work

GE, imagination at work, and GE monogram are trademarks of General Electric Company.

ÄKTA, HiTrap, Sepharose, Rainbow and UNICORN are trademarks of GE Healthcare companies.

Cibacron is a trademark of Ciba-Geigy Corp.

Any use of UNICORN software is subject to GE Healthcare Standard Software End-User License Agreement for Life Sciences Software Products. A copy of this Standard Software End-User License Agreement is available on request.

© 2013 General Electric Company – All rights reserved. First published Dec. 2013.

All goods and services are sold subject to the terms and conditions of sole of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.

GE Healthcare UK Limited Amersham Place Little Cholfont Buckinghamshire, HP7 9NA UK GE Healthcare Europe, GmbH Munzinger Strasse 5 D-79111 Freiburg Germany GE Healthcare Bio-Sciences Corp. 800 Centennial Avenue, P.O. Box 1327 Piscataway, NJ 08855-1327 USA GE Healthcare Japan Corporation Sanken Bldg., 3-25-1, Hyakunincho Shinjuku-ku, Tokyo 169-0073 Japan