Porcine Serum, New Zealand Origin

HyClone[™] Porcine Serum is sourced from New Zealand and filtered in the United States. This product is sterile filtered through multiple 200 nm (0.2 µm) pore size-rated filters and is gamma irradiated at 25 to 40 kGy after filtration. Before dispensing, each lot of serum is pooled using true pool technology to ensure uniformity and consistency between bottles (Fig 1).

Key features of Porcine Serum, New Zealand Origin:

- Complete traceability back to original source
- Virus panel testing according to 9 CFR 113.53
- Produced using true pool technology for uniformity and consistency from bottle to bottle

Product storage and handling

Serum should be stored at -10°C or lower. Once thawed, serum should be stored at 2°C to 8°C for up to six weeks in order to maintain quality. If the serum needs to be stored longer than six weeks after opening, it is recommended to aliquot the serum into convenient volumes and refreeze. Handle bottles that have been stored in freezer carefully. Avoid large temperature shifts and protect the serum from exposure to light. Refer to safety data sheet for any safety recommendations. Storage requirements are listed on the product label.

Thawing

Remove serum from storage at -10°C or lower and place in a refrigerator overnight at 2°C to 8°C. Transfer the serum to a 37°C water bath, agitate periodically to mix the solutes concentrated at the bottom of the container. Do not hold the serum at 37°C any longer than necessary after thawing. Thawing serum in a bath above 40°C without mixing can denature the concentrated proteins in the bottom of the container and precipitates might form in the bottle. Thawing serum at higher temperatures is not recommended.

Alternatively, bottles can be placed directly from storage at -10°C or lower into a 37°C water bath. Bottles should be agitated to



Fig 1. Porcine Serum, New Zealand Origin.

enhance mixing and thawing. Turbidity and flocculent material might be present after thawing or after prolonged storage.

Experience indicates that regardless of the method used to thaw serum, it is critical that it is mixed during the thawing process to prevent the formation of gradients and subsequent precipitation. Because of differences in thawing rates of different components, serum will form a gradient if it is not mixed as it thaws. If serum is allowed to remain in such a gradient state, precipitation is likely to occur.

General culture recommendations

Supplementation of classical media such as Dulbecco's Modified Eagle's Medium (DMEM) is recommended at a range between 5% and 10% Porcine Serum, New Zealand Origin to support culture of a wide variety of cell lines and applications.



Quality control testing

Porcine Serum, New Zealand Origin is assayed for gamma globulin, alkaline phosphatase, lactate dehydrogenase, glutamic pyruvic transaminase (SGPT), glutamic oxaloacetic transminase (SGOT), pH, total protein, albumin, blood urea nitrogen, creatinine, total bilirubin, sodium, potassium, calcium, chloride, inorganic phosphorus, osmolality, and glucose. Assays are subject to change without notice.

Test specifications are listed in Table 1. Results are provided as a courtesy for information only.

Table 1. Test specifications

Endotoxin (Limulus amebocyte lysate gel clot assay)	≤ 100 EU/mL
Hemoglobin (spectrophotometric)	≤ 50 mg/dL
Sterility testing (current USP) Bacteria and fungi	No growth
Virus testing (9 CFR 113.53) Fluorescent antibody Transmissable gastroenteritis virus Porcine adenovirus Porcine parvovirus Bovine viral diarrhea virus Rabies Reovirus Cytopathogenic agents (e.g., IBR) Hemadsorbing agents (e.g., PI3)	Not detected Not detected Not detected Not detected Not detected Not detected Not detected
Mycoplasma Large volume, direct culture Hoechst™ DNA stain	Not detected Not detected
Ouchterlony (double immunodiffusion)	Positive for porcine Negative for equine, bovine, and human

Related products HyClone classical media

HyClone classical media are manufactured using ISO 9001- and ISO 13485-certified processes. All raw material components have passed strict quality control testing to ensure the appropriate level of quality. The classical media are hydrated using purified process water and have undergone 0.1 μ m sterile filtration.

HyClone phosphate buffered saline (PBS)

Our PBS products are manufactured using ISO 9001- and ISO 13485-certified processes. The products have full traceability and documented origin of all formula ingredients.

HyClone trypsin protease

Our trypsin protease is derived from porcine pancreas and is gamma irradiated prior to hydration and filling. The product is formulated without calcium and magnesium.

Ordering information

Size	Product code
500 mL	SH30908.03
1000 mL	SH30908.04
	500 mL

Find certificates of analysis, safety data sheets, standard formulations, product inserts, and protocols at cytiva.com/hyclonecerts.

cytiva.com/hyclone

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate. HyClone is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

Hoechst is a trademark of Hoechst GmbH. All other third-party trademarks are the property of their respective owners.

© 2020 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

