

Biologic manufacturing capacity expansion with single-use technologies

Key variables to consider



The following presentation is based on a simulation of a 2×2000 L mAb process. The simulation compares a single-use process train (SU) with a comparable stainless steel-based process train (SS), both modeled in a traditional, stick-built facility. Are you interested in a simulation of your biomanufacturing process? With our range of simulation tools, we can assist in this, whether it is for scale up/down, "de-bottlenecking", process intensification, transition from stainless to single use, or general optimization.



Manufacturing setup simulation
for 2 × 2000 L mAb process

Titer	3 g/L
Overall process yield	70%
Number of products per year	4
Facility utilization	80%
mAb output per year	168 kg

	TITER		
	2.0	3.0	4.0
2 × 500 L	28	42	56
2 × 1000 L	56	84	112
2 × 2000 L	112	168	224
4 × 2000 L	224	336	448


Assumptions: 70% recovery through purification of 20 batches per year, per reactor




Time to market and capital expenditure

Single-use vs stainless steel technologies

Time to market 

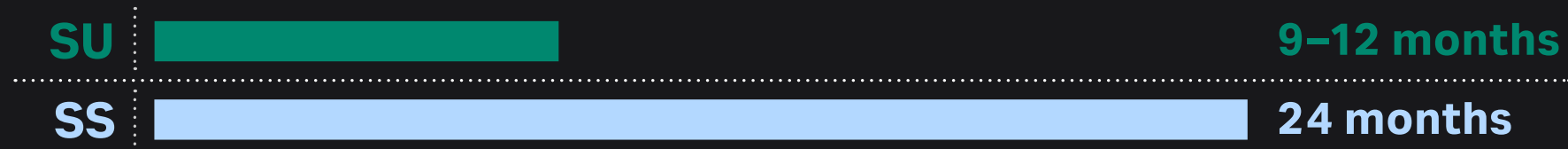

CAPEX 

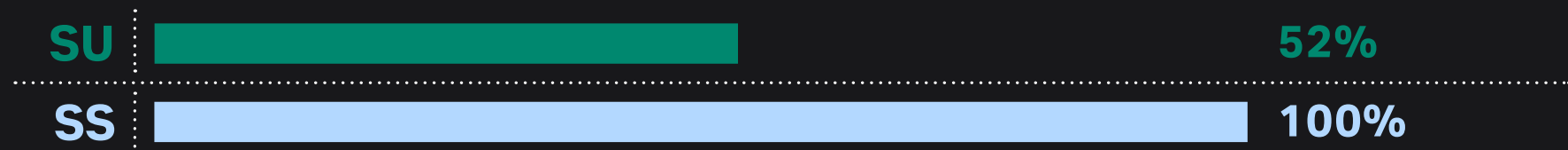

CAPEX = capital expenditure



Stainless steel technologies
take longer to procure, source, qualify, and validate.
The initial cost is also higher.




Single-use technologies
provide faster time to market at lower capital expenditure.


Operating expense

Single-use vs stainless steel technologies

OPEX 

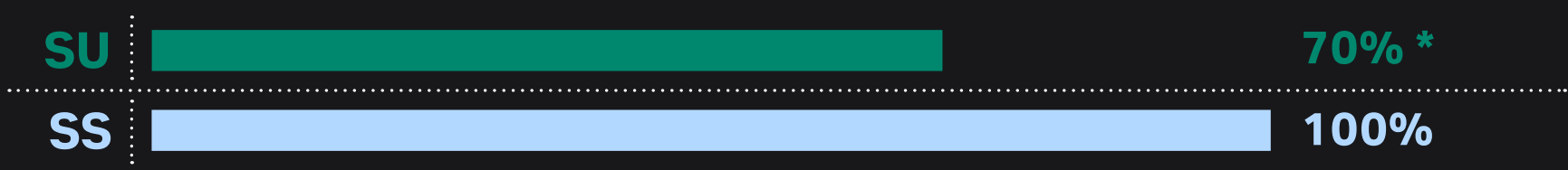

OPEX — labor 

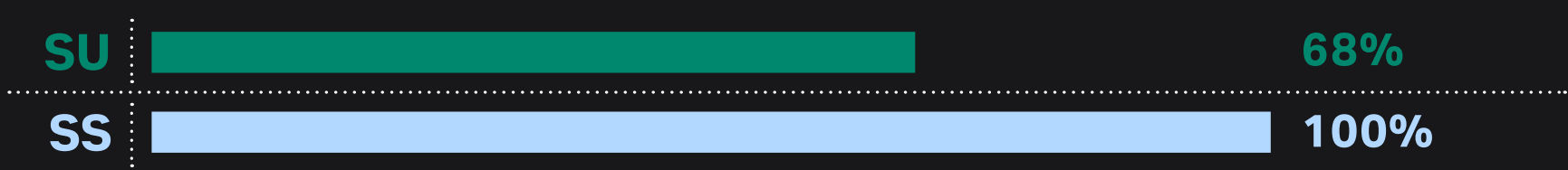

* Industry reports vary from 70% to 120%

OPEX = operational expenditure




Single-use technologies allowed 32% reduction in labor, based on elimination of cleaning and sanitization in place (CIP, SIP), and related testing.



Changeover and output

Single-use vs stainless steel technologies

Changeover 

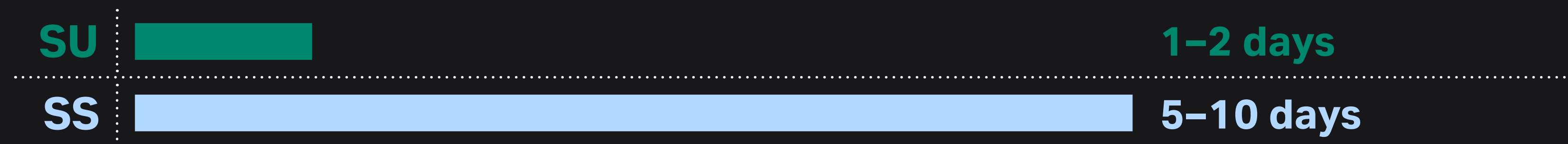


Output 

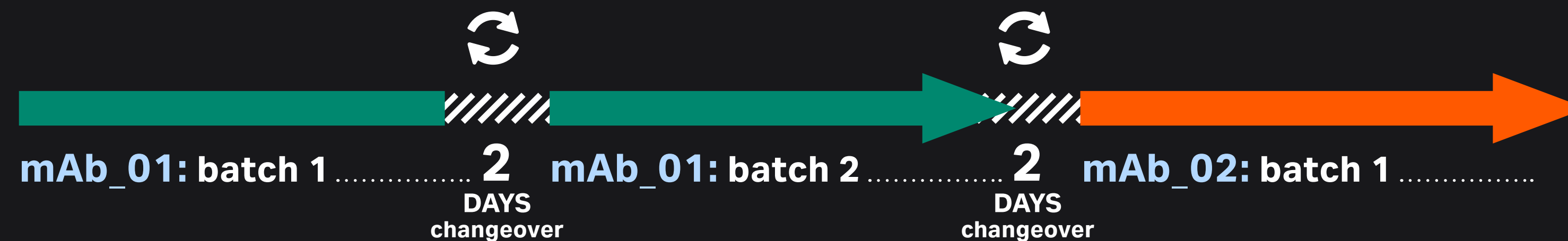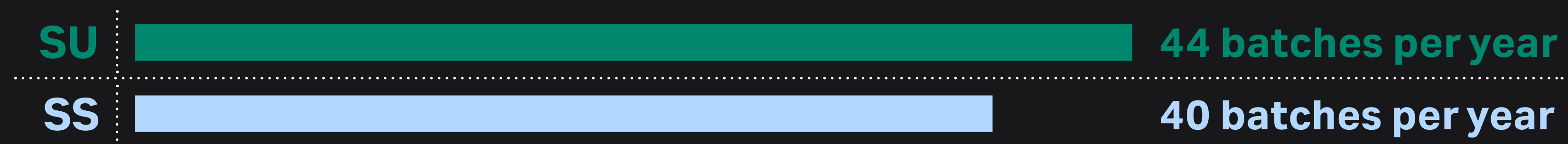



Footprint

Single-use vs stainless steel technologies

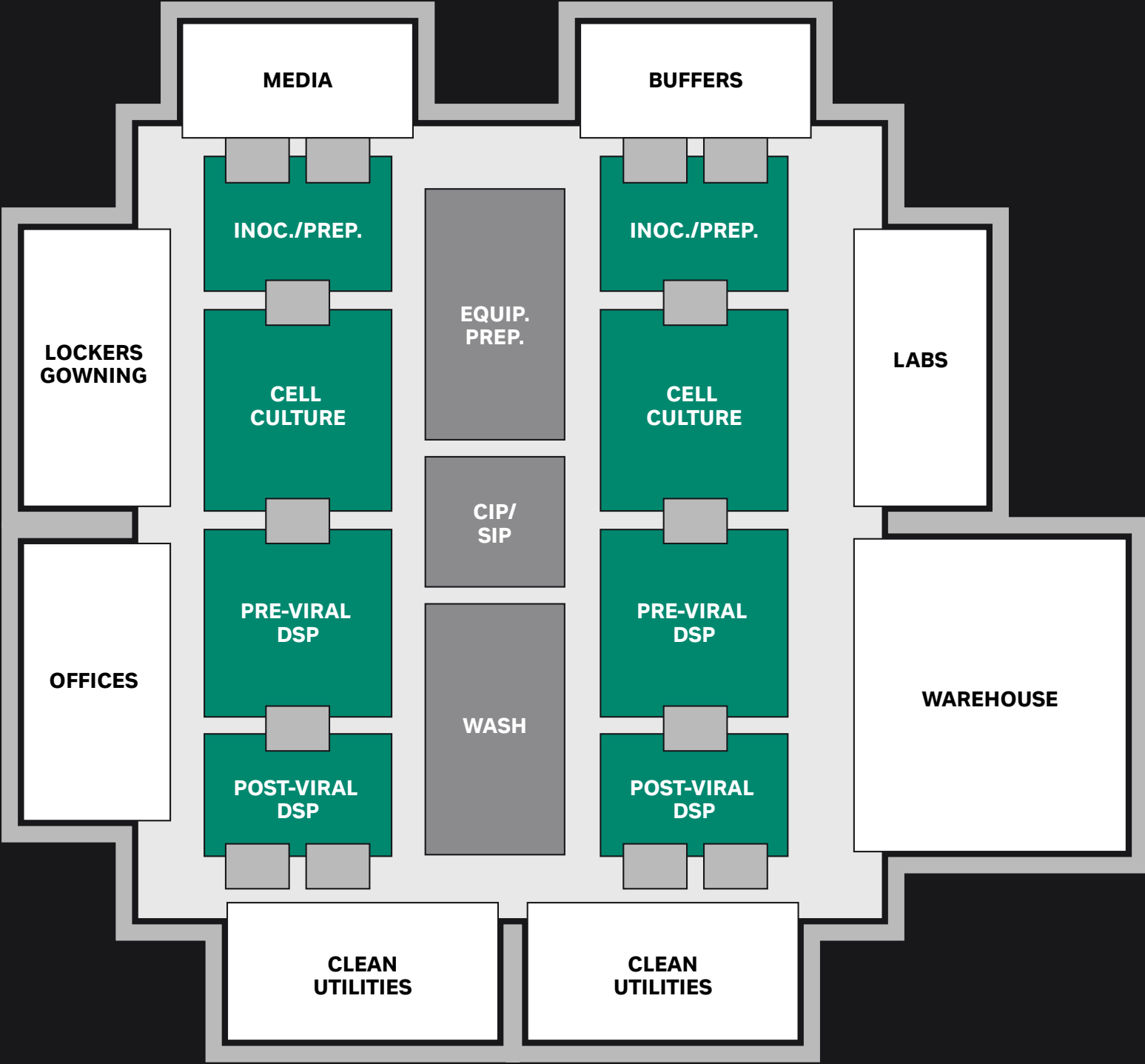

Stainless steel

Footprint 

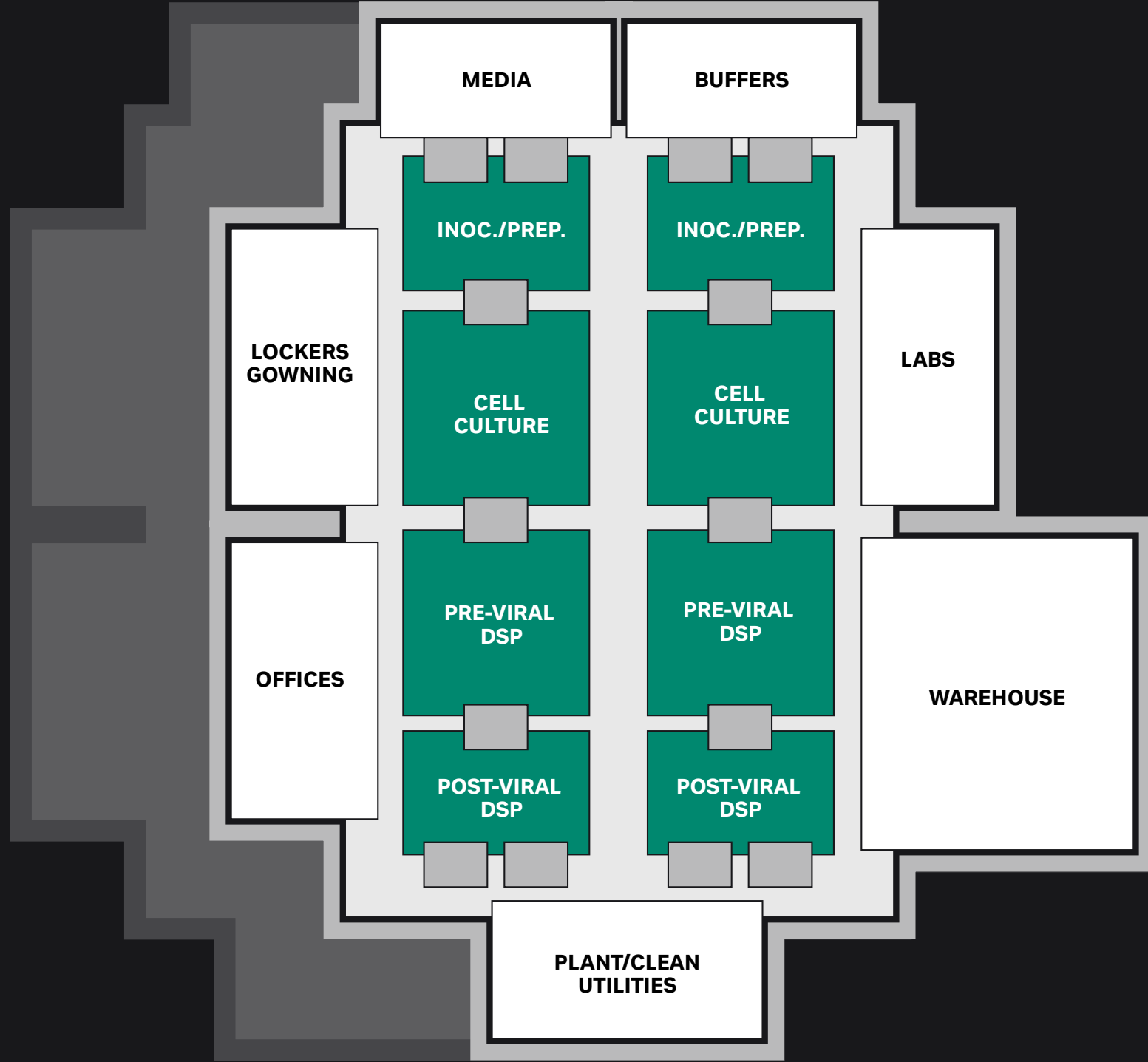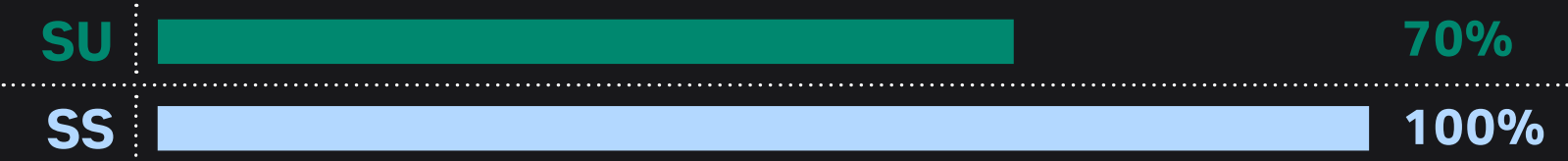


Single-use

Single-use technologies enable smaller facility, less cleanroom space — resulting in lower utilities and HVAC costs.

Cytiva's single-use technologies

Across the entire bioprocess workflow

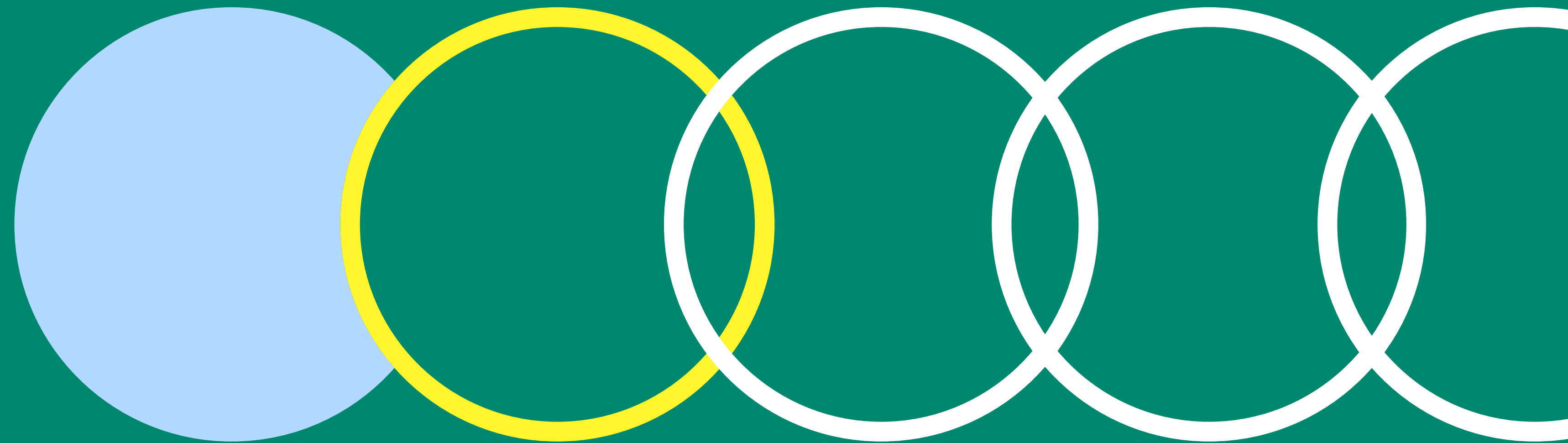


Medium preparation:

Xcellerex™ XDUO 100 to 2500 L mixers, HyClone™ cell culture media

Buffer preparation:

Xcellerex XDUO 100 to 2500 L mixers, HyClone buffers and process liquids

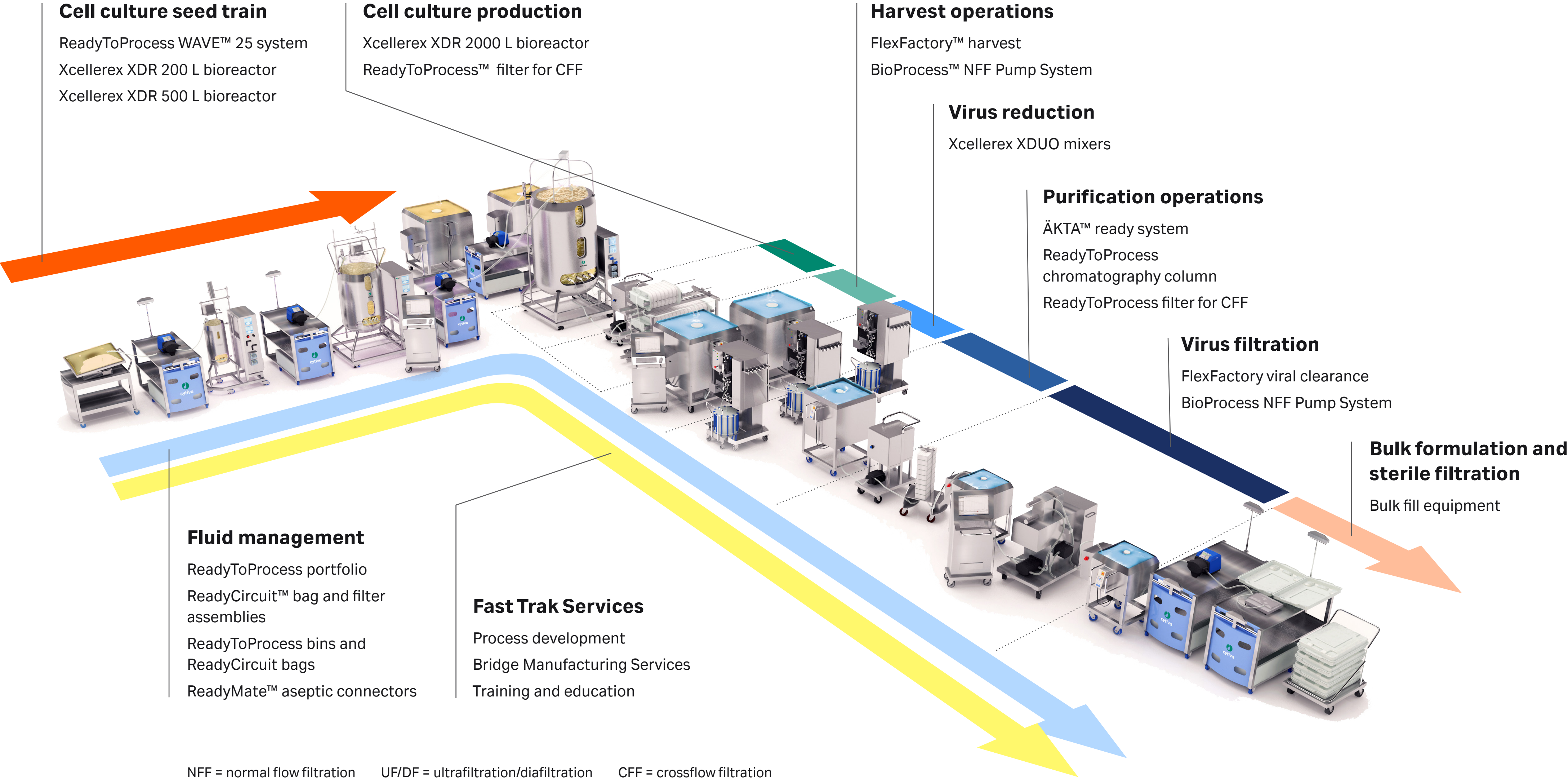

cytiva.com/singleuse

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate. ÄKTA, BioProcess, FlexFactory, HyClone, ReadyCircuit, ReadyMate, ReadyToProcess, ReadyToProcess WAVE, and Xcellerex are trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

© 2020 Cytiva
For local office contact information, visit cytiva.com/contact

CY14204-11Sep20-BR

