

Selection guide

Maximize performance with HyClone media and supplements



Xcellerex XDR-1000

Boost performance Strengthen results

Intensify your cell culture process

Achieving peak performance of protein-producing cell lines in biopharmaceutical manufacturing requires careful selection of cell culture media. Medium supplements and optimally designed feeding strategies can further improve titers and deliver desired product quality characteristics. To help you maximize cell culture performance in your biomanufacturing process, we offer a comprehensive range of HyClone™ media and feeds for many different industrial cell lines, including Chinese hamster ovary (CHO) and human embryonic kidney 293 (HEK293) cells. Our products take you from transfection, screening, and development, to large-scale manufacturing.

Off-the-shelf solutions for cell culture

Media for monoclonal antibody (mAb) and recombinant protein production

These serum-free basal media are animal-derived component-free (ADCF), chemically defined (CD), and protein-free formulations designed to be used with common protein-producing cell lines, such as CHO and HEK293 cells (Table 1).

Table 1. Composition of serum-free media (SFM) for mAb and recombinant protein production

Cell type	SFM	ADCF	CD	Protein-free [‡]	Recombinant protein	Growth factors (peptides)	Hydrolysates	Hypoxanthine/thymidine	Lipids/cholesterol	Poloxamer 188	Product code powder	Product code liquid
CHO*	ActiPro™	•	•	•							Powder and liquid	SH31037 SH31039
	ActiSM™	•	•	•							Powder and liquid	SH31038 SH31040
	CDM4CHO	•	•	•		•			•		Liquid	SH30556 SH30557/SH30558
	HyCell™ CHO	•	•			•		•	•		Liquid	SH30933 SH30934
	SFM4CHO				•	•	•		•		Liquid	SH30518 SH30549/SH30548
	SFM4CHO-Utility					•	•		•		Liquid	SH30517 SH30516
	PF-CHO LS				•	•	•		•		Liquid	NA SH30359
	PF-CHO MPS				•	•	•		•		NA	SH30333 NA
	CDM4NS0	•	•	•					•	•	Liquid	SH30578 SH30579
	CDM4PERMAb	•	•	•		•			•	•	Liquid	SH30872 SH30871
HEK293	CDM4HEK293	•	•	•		•					Liquid	SH30859 SH30858
Hybridoma/myeloma[†]	ADCF-MAb	•			•		•	•	•		Liquid	SH30635 SH30349/SH30547
	CDM4MAb	•	•		•			•	•		Liquid	SH30800 SH30801/SH30802
	CDM4NS0	•	•	•				•	•		Liquid	SH30578 SH30579
	SFM4MAb						•	•	•		Powder and liquid	SH30535 SH30391/SH30513
	PF-MAb	•	•	•							NA	NA SH30138
PER.C6™	CDM4PERMAb	•	•	•		•		•	•		Liquid	SH30872 SH30871

* CHO cell media are for CHO-K1, CHO-M, CHO-S, DG44, DUXB11, GS-CHO and other CHO-derived cell lines

[†] Hybridoma/myeloma media is for cell lines such as NS0, Sp2/0 and P3-derived hybridomas

[‡] Protein-free media do not contain any proteins of molecular weight > M_r 10 000.

Medium supplements

HyClone Cell Boost™ supplements are used for feeding in recombinant protein production to enhance product titer and protein quality. The supplements are designed to provide nutrient formulations that meet your cell line's specific requirements (Table 2).

Basal media and Cell Boost combination

Cell Boost 1 to 7b were screened using a DoE-based approach to select the best performing combination in batch cultivations using various HyClone basal media and CHO cell lines. The general recommendations on basal media and Cell Boost supplements should constitute a starting point for further optimization of feed regimes. The study showed that a combination of Cell Boost 1, 2, 3, 4, 7a, and 7b are likely to fit a broad range of CHO cell lines (Table 3). More information on this study can be found in the application note Optimization of fed-batch culture conditions for a mAb-producing CHO cell line (KA4131090718AN).

Table 2. Composition of Cell Boost supplements

Supplement	Cell type	Amino acids	Vitamins	Glucose	Trace elements	Growth factors (peptides)	Hypoxanthine/thymidine	ADCFLipids	ADCF cholesterol	Product code
Cell Boost 1	CHO, HEK293	•	•	•						SH30584
Cell Boost 2	CHO, PER.C6	•		•						SH30596
Cell Boost 3	Hybridoma, myeloma	•	•	•	•		•			SH30825
Cell Boost 4	CHO	•	•	•	•	•		•	•	SH30857
Cell Boost 5	CHO, HEK293, Hybridoma, NS0	•	•	•	•	•	•	•	•	SH30865
Cell Boost 6	CHO, HEK293, Hybridoma, NS0, T-cells	•	•	•	•	•	•	•	•	SH30866
Cell Boost 7a	CHO	•	•	•	•					SH31026
Cell Boost 7b	CHO	•								SH31027

Table 3. The suitability for use of selected Cell Boost supplements with various CHO cell lines

	Basal media	Cell Boost 1	Cell Boost 2	Cell Boost 3	Cell Boost 4	Cell Boost 7a	Cell Boost 7b
CHO-S (mAb 7)	ActiPro*					+	+
	CDM4PERMAb	+	+		+	+	+
	CDM4MAb	+	+		+		
DG44 (mAb 5)	ActiPro	+				+	+
	CDM4NSO	+	+	+		+	+
CHO-M	ActiPro	+		+		+	+
	CDM4NSO	+		+		+	+

+ Good performance

* No screening of Cell Boost 1 to 7b performed: recommendation on Cell Boost 7a and 7b for ActiPro is based on product description and previous studies.

Table 4. Composition of media for vaccines and viral vectors

Cell type	SFM	ADCF	CD	Protein-free [‡]	Recombinant protein	Growth factors (peptides)	Hydrolysates	Hypoxanthine/thymidine	Lipids/cholesterol	Poloxamer 188	Product code powder	Product code liquid	
EB66	CDM4Avian	•	•	•	•						Liquid	SH31035	SH31036
HEK293	CDM4HEK293	•	•	•		•					Liquid	SH30859	SH30858
	SFM4HEK293	•		•	•		•		•		Liquid	SH30522	SH30521
	PF-293	•		•		•	•		•		Liquid	SH30355	SH30356
Sf9, Sf21, High Five™	SFM4Insect	•					•		•		Powder and liquid	SH30912	SH30913
	SFX-Insect						•		•		Liquid	SH30350	SH30278
	TMN-FH						•				NA	NA	SH30280
PER.C6	CDM4PERMAb	•	•	•		•		•	•		Liquid	SH30872	SH30871
	CDM4Retino	•	•	•		•		•	•		Liquid	SH30519	SH30520
Vero, COS-7, MDCK, MDBK	SFM4MegaVir	•		•		•	•		•		NA	SH30587	NA

[‡] Protein-free media does not contain any proteins of molecular weight > M_r 10 000.

Table 5. Composition of transfection media

Cell type	SFM	ADCF	Recombinant protein	Growth factors (peptides)	Hypoxanthine/thymidine	Lipids/cholesterol	Poloxamer 188	Product code powder	Product code liquid
CHO	HyCell TransFx-C [†]	•	•	•	•	•	NA	SH30942	SH30941
HEK293	HyCell TransFx-H [†]	•	•	•	•	•	NA	SH30944	SH30939
	SFM4Transfx-293 [‡]	•	•	•			Liquid	SH30861	SH30860

[†] Developed for transient transfection and recombinant protein production.

[‡] Developed for stable transfection and transfection of lentiviral and adenoviral constructs and production of virus.

Media for vaccine and viral vector production

Our portfolio includes media for Vero, MDCK, MDBK, and COS-7 cells, for production of vaccines against, for example, influenza, polio, and MMR, as well as for EB66® and other cell lines used in the production of viral vectors (Table 4).

Media for transfection and transient expression

Our transfection media have been tested with a wide range of HEK293 and CHO cell lines and support high transfection efficiency using lipid-mediated (e.g., DharmaFECT™), polymer-mediated, and other transfection methods (Table 5). HEK293 transfection medium is suitable for adenovirus (AdV), adeno-associated virus (AAV), lentivirus, retrovirus, and recombinant protein production. CHO cell transfection medium is suitable for recombinant protein and mAb production.

About our media and supplements for bioprocessing applications

For your convenience, our media are available in both liquid and powder format in a variety of package sizes, which can be customized to meet your needs. To support your large-scale manufacturing needs, the media can be custom manufactured in lot sizes up to 10 000 L for liquid and 6500 kg for powder (density-dependent). If you prefer a customized medium formulation, ask about our Fast Trak medium development and optimization services.

Cytiva is a global provider of biomanufacturing solutions. Our medium and supplement manufacturing operations are part of our holistic security of supply program, based on the three pillars of supply chain sustainability, business continuity, and communication.

cytiva.com

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate. ActiPro, ActiSM, Cell Boost, DarmaFECT, HyCell, and HyClone are trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

EB66 is a registered trademark of Valneva. High Five is a trademark of Thermo Fisher Scientific. PER.C6 is a trademark of Janssen Vaccines and Prevention B.V. All other third-party trademarks are the property of their respective owners.

© 2020 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit [cytiva.com/contact](https://www.cytiva.com/contact)

CY16150-10Dec20-SG

