

# CPAK-100 and CPAK-101 Sampling Lines

## CELL PROCESSING


The sampling lines CPAK-100 and CPAK-101 are single-use accessories for retrieving product sample. With a sterile connecting device (SCD) connect a sampling line to your product bag or line of any single-use kit; then use CPAK-100 for samples up to 3 mL and CPAK-101 for samples up to 8 mL. The sampling lines are supplied sterile and are equipped with a microbial filter ensuring their sterility when properly used.

### Key features

Part	Name	Description
1	Tubing compatible with SCD connection	PVC tubing with an external diameter of 4 mm, compatible with SCD connection
2	Sampling pillow	Graduated pillow with capacity up to 3 mL for CPAK-100 and up to 8 mL for CPAK-101
3	Needleless port	Check valve with cap to connect a syringe with luer port

### Product specifications

Shelf life	Two years from the date of manufacture
Sterility	Ethylene oxide, non-pyrogenic
Material	Standard medical grade PVC, PTFE, silicone and PP

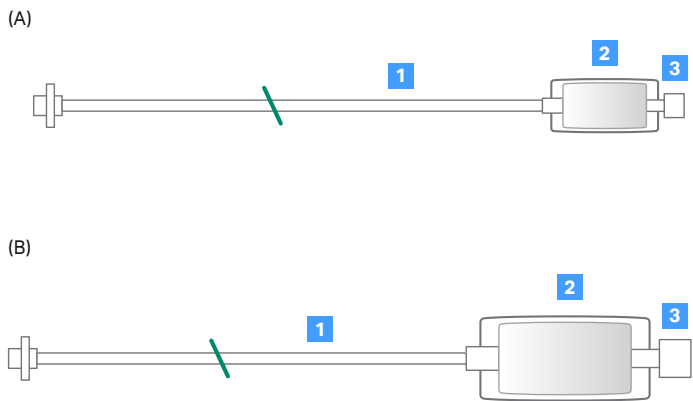

**Fig 1.** (A) CPAK-100 Sampling Line with a capacity of 3 mL, (B) CPAK-101 Sampling Line with a capacity of 8 mL.

### Storage requirements

Store sampling lines in a clean and dry environment without chemical or biological contamination within the following conditions:

	Operation	Storage and transport
Temperature	7°C to 27°C	4°C to 40°C
Relative humidity	30% to 75%, non-condensing	20% to 75%, non-condensing

# Ordering information

Product	Product code
CPAK-100 Sampling Line	29275109
CPAK-101 Sampling Line	29275108
<ul style="list-style-type: none"><li>• Each unit is packed individually</li><li>• Minimum order: Box of 60 units (dimensions: 33 × 35 × 13 cm, 13 × 14 × 5 in)</li></ul>	

Related products	Product code
Sepax™ C-Pro Cell Processing instrument	29264741
CT-49.1 Sepax C-Pro Cell Processing kit	29264738
CT-60.1 Sepax C-Pro Cell Processing kit	29264739
CT-90.1 Sepax C-Pro Cell Processing kit	29264740
Sefia™ S-2000 Cell Processing Instrument	29285527
CT-300.1 Sefia Cell Processing Kit	29284866
CT-800.1 Sefia Cell Processing Kit	20001

## cytiva.com/celltherapy

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate. Sefia and Sepax are trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

© 2020 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit [cytiva.com/contact](https://cytiva.com/contact)

CY14614-18Sep20-DF

