

mRNA manufacturing workflow

Plasmid linearization



Objective: Produce DNA template

Considerations

- Restriction enzyme digest of plasmid DNA (pDNA) to create a linear template
- Quality of ingoing pDNA to control isomeric form

Strategies

- Choose appropriate scale and vessel with controlled environment
- Select plasmid source and use analytics with quality control in mind

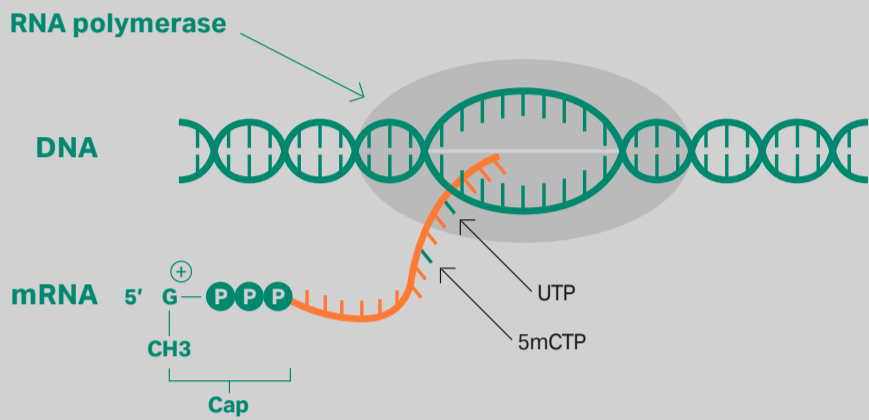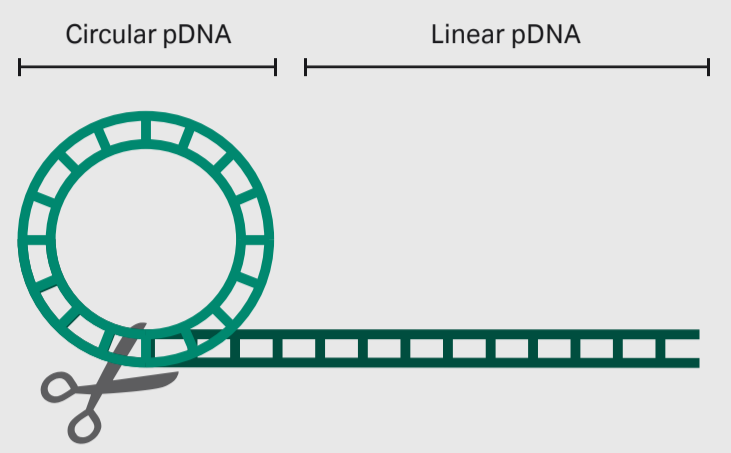

In vitro transcription (IVT)

Objective: Enzymatic synthesis and capping of mRNA

Considerations

- Choice of enzymes, dNTPs and nucleosides
- Whether to cap before or during transcription
- DNase treatment and removal of residual DNA and RNA isoforms
- Scale and use of organic solvents

Strategies

- Identify enzymes and nucleotides specific to target mRNA; replace uridine and/or cytosine with modified nucleosides
- Precipitation and wash step to remove residual impurities
- Optimize reaction conditions for capping (e.g., temperature)

Buffer exchange

Objective: Prepare feed for purification step


Considerations

- Avoiding loss of titer for best product recovery
- Mitigating risk of failure in multistep process
- Avoiding filter fouling, clogging, and contamination


Strategies

- Select the right filter type and sizing for scale-up
- Use a scalable, automated TFF system
- Consider single-use, closed systems

ÄKTA readyflux™ system with ReadyToProcess™ hollow fiber cartridges

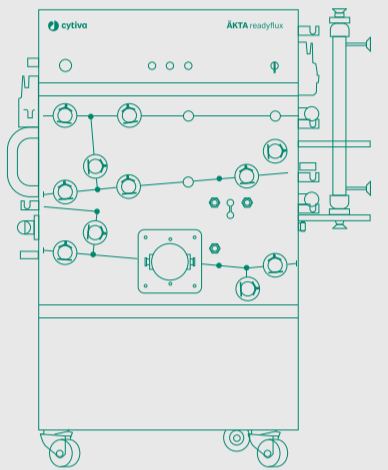

Oligo(dT) resins

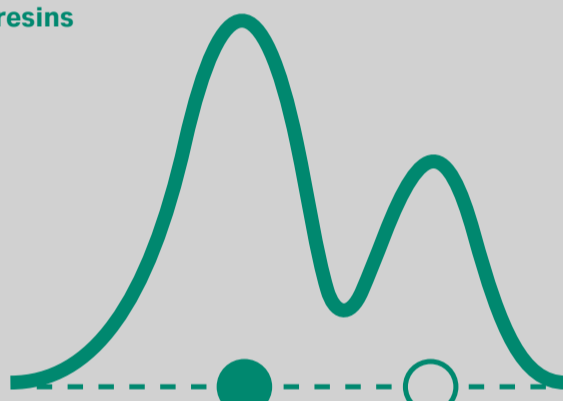

Capture

Objective: mRNA purification

Considerations

- Molecule size – mRNA is ~10x larger than proteins
- Selecting resin that purifies full-length mRNA + removes contaminants
- Elution conditions for maximum recovery

Strategies

- Use resins with specific ligand for faster throughput (promising option: fiber-based chromatography)
- Control bioburden + speed changeover with prepacked resin formats
- Use scalable column platform from PD to full-scale manufacturing

Concentration

Objective: Volume reduction and change to stable salt conditions


Considerations

- Avoiding loss of titer for best product recovery

Strategies

- Choose a scalable, automated, single-use, closed TFF system
- Tailor buffer conditions to control secondary mRNA structures
- Optimize concentration with filter cut-off

ÄKTA readyflux™ system with ReadyToProcess™ hollow fiber cartridges

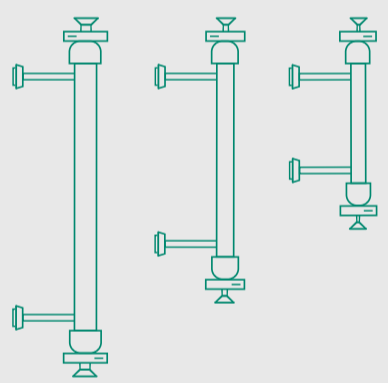

LNP formation

Objective: Lipid nanoparticle (LNP) formation as an mRNA drug delivery vehicle

Considerations

- Optimizing mix of mRNA with lipids to generate desired LNP size
- Optimizing ratio of empty vs filled LNPs
- Avoiding excess lipid monomer formation

Strategies

- Monitor particle size, polydispersity index, LNP formation efficiency
- Validate chemistry, molecular weight cut-off/pore size of filters
- Validate parameters (e.g., flow rate, path length, shear rate, shear stress)
- Ensure scalability – optimized parameters from bench to process scale

Polishing

Objective: Remove product and process impurities


Considerations

- Optimal balance – recovery of encapsulated mRNA and removal of free lipids and non encapsulated mRNA
- Matching polishing solution to impurity type

Strategies

- Choose core beads to scavenge impurities and meet regulatory demand
- Tangential flow filtration with appropriate cut off for optimal recovery and efficient impurity removal

**ReadyToProcess™ hollow fiber cartridges
Capto™ Core 700 chromatography resin
ÄKTA ready™ single-use chromatography system**

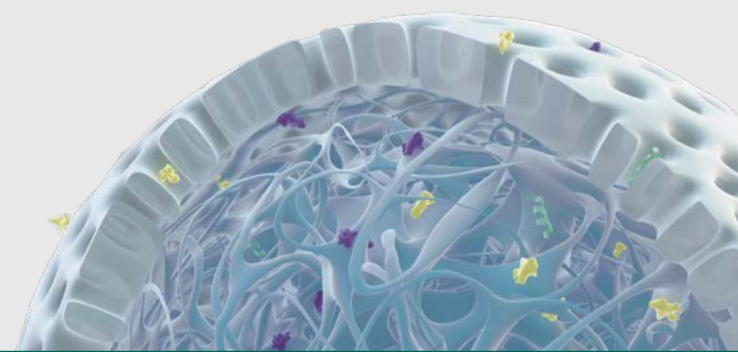

Drug product

Objective: Drug product for delivery to patient

Considerations

- Which path? Personalized or platform mRNA drug product
- Long-term storage conditions
- Visual inspection of translucent drug product

Strategies

- Scale out to add capacity quickly, especially in personalized
- Co-locate drug product with drug substance to streamline production and reduce supply chain risk
- Options to transition from vial to syringe for launch

SA25 Aseptic Filling Workcell



[cytiva.com/mrna](https://www.cytiva.com/mrna)

Cytiva and the Drop logo are trademarks of Life Sciences IP Holdings Corp. or an affiliate doing business as Cytiva. ÄKTA ready, ÄKTA readyflux, Capto, and ReadyToProcess are trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

© 2021 Cytiva
CY19802-26Oct21-IG
GL-21-00354

Learn more about mRNA manufacturing [here](#).