



# ImageQuant™ TL 11.0

## License Setup Guide for macOS

# Table of Contents

- 1 Introduction ..... 3**
- 2 Collect license using activation ID ..... 4**
- 3 Administer electronic license ..... 7**
- 4 Product support ..... 8**

# 1 Introduction

## About this document

This document describes the licensing features, license configuration, and management of ImageQuant™ TL (IQTL).

# 2 Collect license using activation ID

## Introduction


The ImageQuant TL software requires an electronic license (eLicense) to operate. The license file is collected from the Cytiva software licensing web page using an activation ID for Mac computers.

The activation ID is delivered via e-mail after ordering ImageQuant TL. Store the activation ID in a safe place.


## Collect and activate the electronic license

Follow the steps below to collect and activate the electronic license.

- | Step | Action   |
|------|--|
| 1    | Go to <a href="https://cytiva.com/eDelivery">cytiva.com/eDelivery</a> and log in by selecting <b>Activation ID</b> and typing in the activation ID that you received by e-mail during your purchase. |
| 2    | On the top toolbar, click <b>Activation &amp; Entitlements</b> → <b>List Entitlements</b> .  |

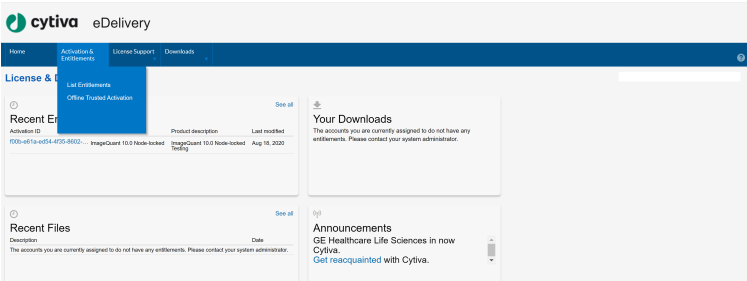

- 3 Select the entitlement in the list for Product name 4, and activate it by clicking **Action** → **Activate**.

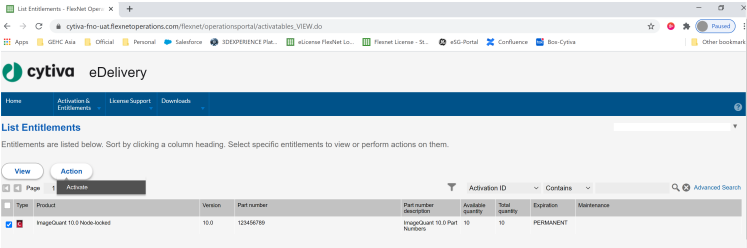


Step      Action

4      Type your personal details and click **Next**.

5      Under **Configure Hosts**, click  to add host.

6      In **Create License Host** window, type the **Host ID** (also called MAC address) for the computer where you have installed ImageQuant TL. Then click **OK** to generate the license file. For details on how to find the Host ID, refer to [Find Host ID \(MAC address\) for macOS, on page 6](#).

Step	Action
7	Under <b>License Details</b> , select <b>View</b> → <b>Save to File</b> to save the license file to any folder you want on your computer.

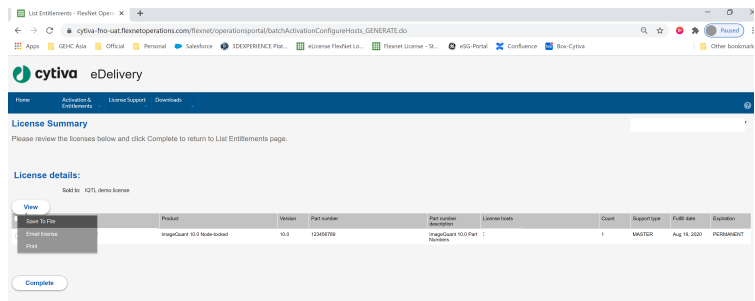

**Result:**  
The license file is ready to be used.

## Find Host ID (MAC address) for macOS

Follow the steps below to find the **Host ID** for macOS.

Step	Action
1	Open the Apple menu and click the Apple icon.
2	Scroll to <b>System Preferences</b> .
3	Select <b>Network</b> → <b>Wi-Fi</b> → <b>Advanced</b> .
4	Copy the MAC address shown under the <b>Wi-Fi Address</b> or <b>Airport Address</b> .

**Note:**  
*If the system only connects via Ethernet, then use the Ethernet address.*

# 3 Administer electronic license

## Introduction


For details on how to collect the eLicense using this activation ID, see [Chapter 2 Collect license using activation ID, on page 4](#). When the eLicense has been downloaded onto your computer, see instructions in the following subsection to activate and begin using the software.

**Note:** *If the activation ID has not been delivered within 7 days of the order, contact technical support.*

## Node-locked licence

Follow the steps below when using a node-locked license.

Step	Action
1	Download and save the node-locked license in any folder on your computer.
2	Start the ImageQuant TL software.
3	When the ImageQuant TL control hub opens, click <b>Configure license</b> .
4	Browse to the folder where you saved the node-locked license and click on the license file and <b>Done</b> .

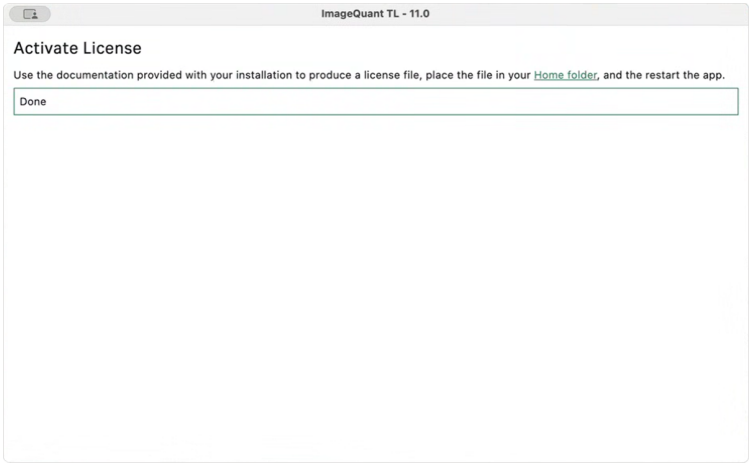

**Result:**  
The software is ready to use.

5	Return to the hub.
---	--------------------

## 4 Product support

### Contact us

If there are any problems with the installation or use of ImageQuant TL, contact your local product support team, or use the contact form found at [cytiva.com/contact](https://cytiva.com/contact).

Page intentionally left blank

**Give feedback on this document**

Visit [cytiva.com/techdocfeedback](https://cytiva.com/techdocfeedback) or scan the QR code.



## cytiva.com

Cytiva and the Drop logo are trademarks of Life Sciences IP Holdings Corporation or an affiliate doing business as Cytiva.

ImageQuant is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

Any other third-party trademarks are the property of their respective owners.

© 2025 Cytiva

ImageQuant © 2025 Cytiva

Any use of software may be subject to one or more end user license agreements, a copy of, or notice of which, are available on request.

For local office contact information, visit [cytiva.com/contact](https://cytiva.com/contact)

29751073 AA V:3 12/2024