

ÄKTA go Unpacking Instructions

Read this before unpacking ÄKTA go



This instruction contains information that is important for safety during the unpacking of ÄKTA™ go. Read the whole instruction carefully before starting to unpack the system.

When you receive the delivery

- Record on the receiving documents if there is any apparent damage to the delivery box. Inform your Cytiva representative of such damage.
- Move the delivery box to a protected location indoors.

About this instruction

This instruction covers the unpacking of ÄKTA go. See *Operating Instructions 29360951* for instructions on how to install the instrument.

The delivery box contains the ÄKTA go instrument and a cardboard box. The cardboard box contains the top and bottom trays, product documentation, and an accessories box.

Instruction

Follow the steps below to unpack the instrument and lift the instrument onto the bench.




CAUTION


Heavy object. Use suitable precautions when moving the system. Two people are recommended to lift the system safely.

Step	Action
------	--------

- | | |
|---|-----------------------------------------------------------------|
| 1 | Cut the plastic straps, and remove the lid and protective foam. |
|---|-----------------------------------------------------------------|

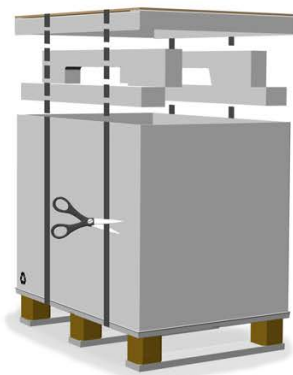


- | | |
|---|---------------------------------------------------------------------------------------------------|
| 2 | Lift off and store the cardboard box that contains ÄKTA go product documentation and accessories. |
|---|---------------------------------------------------------------------------------------------------|

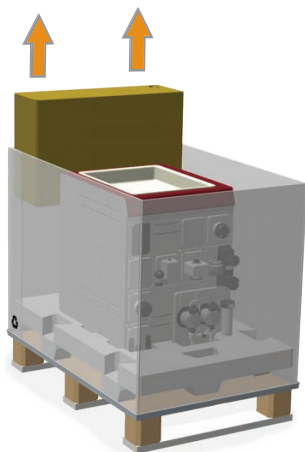

Step	Action
------	--------

- | | |
|---|--------------------------------------------------------------|
| 3 | Remove the cardboard hood and the cover from the instrument. |
|---|--------------------------------------------------------------|

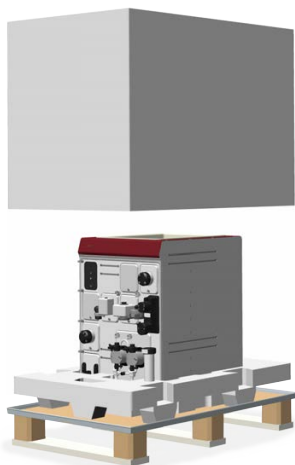

- | | |
|---|---------------------------------------------------------------------------------------------------------|
| 4 | Using two persons, grip the instrument from the sides and carefully lift the instrument onto the bench. |
|---|---------------------------------------------------------------------------------------------------------|

Note:

The instrument is heavier at the front. Do not tip the instrument when lifting.





cytiva.com/AKTA

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

AKTA is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

All other third-party trademarks are the property of their respective owners.

© 2020 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

29383543 AC V:5 12/2020