

## Optimization and Integration of Single-Pass, Continuous Final Formulation Technologies for mAb Purification

Heather MalloBlaine Trafton, Jeannette Marshall; Pall Life Sciences - BioPharm Applications R&D; 20 Walkup Drive, Westborough, MA 01581 USA

### INTRODUCTION

▶ The utilization of Pall's patented Cadence™ single-pass TFF (SPTFF) technologies is a key enabler of continuous bioprocessing initiatives. The integration and optimization of a prototype Cadence Inline Diafiltration (ILDF) module with the Cadence Inline Concentrator (ILC) with sterile filtration is key to development of a fully continuous, streamlined final formulation process platform. Coupling these continuous final formulation processes allows for further optimization and reduction of module/device flushing as well as reduction of in-process buffer consumption compared to independent operation. The implementation of this continuous final formulation train (ILDF/ILC/0.2 μm) successfully achieved > 3-log (> 99.9%) primary buffer species impurity removal and ≥ 20x concentration factor while delivering a 0.2 μm filtered monoclonal antibody.

### OBJECTIVE

- ▶ To provide an operationally simplified data-driven approach and assessment of the SPTFF attributes for the continuous final formulation of mAbs and address the subsequent product and process development requirements.
- ▶ Offer operational and logistical improvements of coupling existing single-pass technologies for maintaining process simplicity and control.

### BACKGROUND

- ▶ Cadence Inline Concentrators are well-suited SPTFF solutions for numerous bioprocess applications and have been shown to remove constraints in existing facilities via:
  - In-process volume reduction applications (i.e., post-harvest, pre-capture, tank-to-tank transfer, etc.)<sup>1-2</sup>,
  - Process coupling runs with affinity, ion-exchange and membrane chromatography steps in batch and continuous mode showcasing enhanced productivity<sup>3-4</sup>,
  - Increased flexibility of manufacturing capabilities with significant cost savings<sup>5-6</sup>, and
  - Achieving high concentration formulations with custom pre-assembled modules available in various flow path configurations<sup>6</sup>.

### RESULTS AND DISCUSSION


#### High Concentration Formulations with Cadence ILC

- ▶ Cadence ILC is an important addition to the process development tool-box for platform process evaluation and is a crucial and proven enabler of integrated, streamlined and continuous bioprocessing initiatives<sup>2-6</sup>.

- ▶ For instance, utilizing ILC for UF2 enables multiple advantages over a conventional UF system, such as:
  - Non-capital expense,
  - Lower system hold-up volumes,
  - High productivity and improved recovery (> 99%), and
  - Reduced over-concentration targets.

**Figure 1**


Evaluation of ILC Volumetric Concentration Factor for Final Formulation Operations (30 kDa T01 module, 0.065 m<sup>2</sup>)

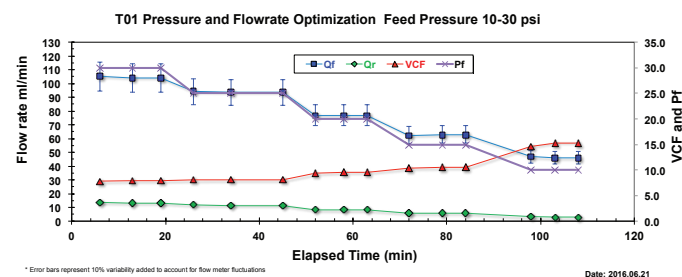


- ▶ Ease of process optimization with ILC enables the generation of a design space achieved with a single module.
- ▶ By assessing the process at a wide range of operating conditions (via feed pressure control) provides a stable and reproducible performance map.

**Figure 2**

Integrated Cadence ILC + 0.2 μm Sterile Filter

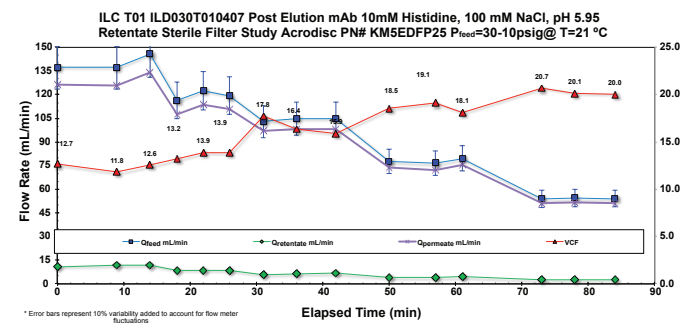

- ▶ Implementation of 0.2 μm sterile filter provides additional back pressure to the Retentate, allowing for increased concentration.
- ▶ Integration provides sterile final product without added equipment and operator interaction for final concentration processes.

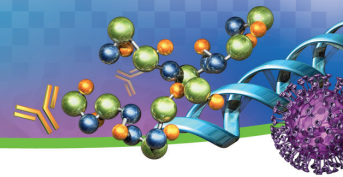


### Continuous, Single-pass and/or Inline Diafiltration (ILDF)

- ▶ ILC design principles and its versatile and robust performance led to the assessment of novel continuous diafiltration strategies, which resulted in a working prototype of a novel inline diafiltration (ILDF) module.
- ▶ Feasibility runs with 10 - 80 g/L hlgG feeds resulted in successful removal efficiencies ( $\geq 99.9\%$  or  $\geq 3\text{-log}$ ) between ILDF process targets and validated test data (Figure 2).


**Figure 3**

Performance benchmark for the continuous, single-pass, inline diafiltration module with 30 kDa regenerated cellulose membrane [Feed: 10 – 80 g/L hlgG in 0.025 M Na acetate + 0.5 M NaCl (47 mS/cm) + 50 g/L Glucose and DF: 0.025 M Na acetate + 0.05 M NaCl (6.7 mS/cm)].

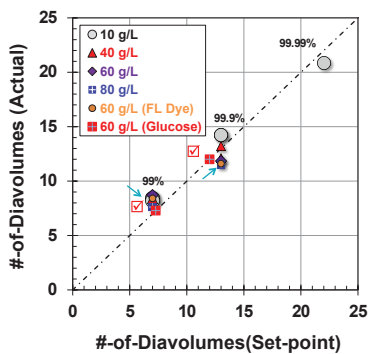

**Figure 4**


Scalability of Cadence ILDF (Feed Pressure vs. LMH)

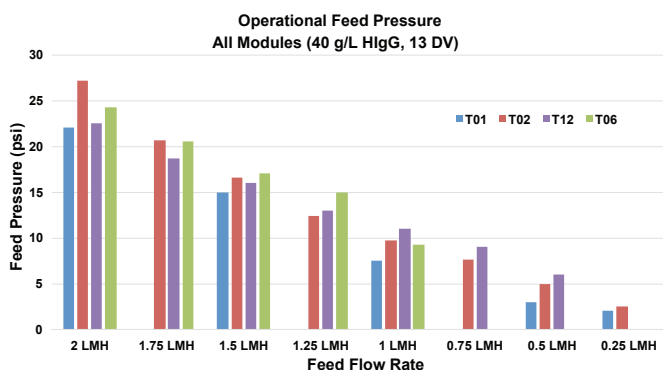

- ▶ ILDF process scalability shown up to >50x scale-up.
- ▶ Scalable feed pressure performance across device flow range providing predictable operational results.


**Figure 5**

Streamlined Continuous Final Formulation Steps with Virus, ILDF, ILC, and Sterile Filtration

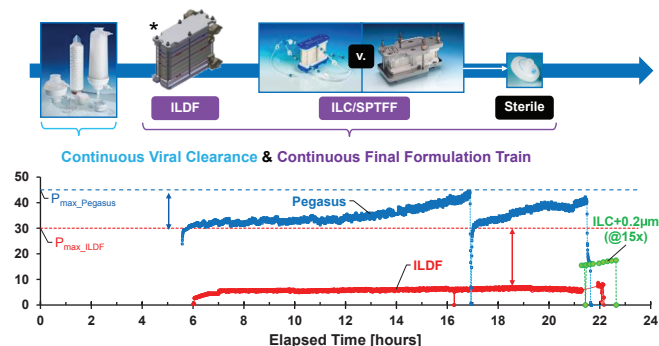

**Figure 6**


Complete Integration of Final Formulation Technologies

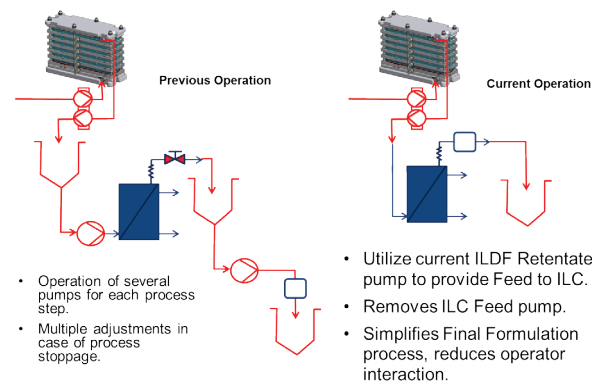


- ▶ Utilization of a single pump to control Feed and Retentate of ILDF maintains concentration throughout processing.
- ▶ ILDF Pump provides feed to the Cadence ILC module without the need for a surge tank or additional pumping equipment.
- ▶ Integrating 0.2  $\mu\text{m}$  sterile filter to the Cadence ILC Retentate removes need for surge tank and additional equipment.
- ▶ Streamlines ease of use operation of SPTFF technologies.

**Figure 7**

Coupled Cadence ILDF-ILC-Sterile Final Formulation Step

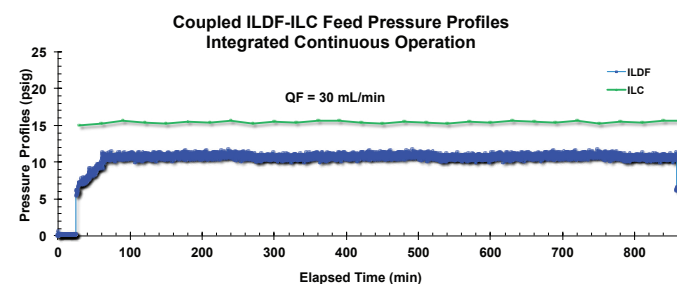

- ▶ Process stability maintained throughout duration of integrated continuous process.

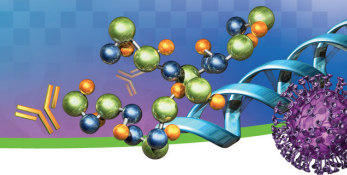

- ▶ Cadence ILC technology platform
  - Provides large operation window and flexibility of concentration factors with minimal control.
  - Presents versatile capabilities for the processing of biologics.
  - Is scalable, flexible and available in various formats and configurations in order to address molecule- and process-specific customization needs.
  - Can remove constraints in existing facilities and increase the flexibility of manufacturing capabilities by increasing productivity and facilitating further use of disposables.
  - Enables process coupling and integrated continuous bioprocessing.
- ▶ The Cadence ILDF design and ease-of process control
  - Addresses an innovation gap for the biopharmaceutical industry.
  - Completes the product/technology offering for the realization of continuous final formulation.
  - Brings the Biopharm industry one step closer to the realization of an end-to-end integrated continuous bioprocessing platform via the integration of ILDF/ILC/ SPTFF technologies.
- ▶ Cadence ILDF-Cadence ILC-Sterile Coupling Integration
  - Provides streamlined approach to final formulation processing.
  - Eliminates multiple surge tanks and operational equipment in between process steps.
  - Reduces operator interaction required for development and process control.

## REFERENCES

- 1 Pall Application Note (USD2789): Cadence™ Systems Employ New Single-Pass TFF Technology to Simplify Processes and Lower Costs (2011).
- 2 Casey, C. and Ayturk, E., Pall Application Note (USTR 2913): Volume Reduction and Process Optimization with Cadence™ Single-Use Inline Concentrator (2013).
- 3 Schofield, M., Rogler, K., Gjoka, X., Ayturk, E., and Gantier, R. Pall Application Note (USD3002): Productivity and Economic Advantages of Coupling Single-pass Tangential Flow Filtration to Multi-Column Chromatography for Continuous Processing (2014).
- 4 Casey, C. Rogler, K., Gjoka, X., Gantier, R. and Ayturk, E. Pall Application Note (USD3003): Cadence™ Single-pass TFF Coupled with Chromatography Steps Enables Continuous Bioprocessing while Reducing Processing Times and Volumes (2015).
- 5 Ayturk, E. and Gantier, R. Process Economics of Continuous mAb Processing with Single-Pass Tangential Flow Filtration. Presented at 249th American Chemical Society National Meeting and Exposition, March 23-26, 2015, Denver, CO, USA.
- 6 Ayturk, E. and Forespring, C. Simplifying Bioprocessing with Single-pass TFF. Genetic Engineering and Biotechnology News, Vol. 36, No. 4, (2016).

### Corporate Headquarters

Port Washington, NY, USA  
+1.800.717.7255 toll free (USA)  
+1.516.484.5400 phone  
biopharm@pall.com e-mail

### European Headquarters

Fribourg, Switzerland  
+41 (0)26 350 53 00 phone  
LifeSciences.EU@pall.com e-mail

### Asia-Pacific Headquarters

Singapore  
+65 6389 6500 phone  
sgcustomerservice@pall.com e-mail

---

Visit us on the Web at [www.pall.com/biopharm](http://www.pall.com/biopharm)


E-mail us at [biopharm@pall.com](mailto:biopharm@pall.com)

---

### International Offices

Pall Corporation has offices and plants throughout the world in locations such as: Argentina, Australia, Austria, Belgium, Brazil, Canada, China, France, Germany, India, Indonesia, Ireland, Italy, Japan, Korea, Malaysia, Mexico, the Netherlands, New Zealand, Norway, Poland, Puerto Rico, Russia, Singapore, South Africa, Spain, Sweden, Switzerland, Taiwan, Thailand, the United Kingdom, the United States, and Venezuela. Distributors in all major industrial areas of the world. To locate the Pall office or distributor nearest you, visit [www.pall.com/contact](http://www.pall.com/contact).

The information provided in this literature was reviewed for accuracy at the time of publication. Product data may be subject to change without notice. For current information consult your local Pall distributor or contact Pall directly.

© 2017 Pall Corporation. Pall, , and Cadence are trademarks of Pall Corporation are trademarks of Pall Corporation. ® indicates a trademark registered in the USA and TM indicates a common law trademark.